Chest freezer ULT C300

220-240V/50Hz

- Low cost of ownership
- Limited maintenance and on-site service
- Easily installable & movable

- Efficient cyclopentane insulation
- Eco-friendly, HCFC-free Nature R refrigerant
- ISO ETL, ATEX Certifications

Specifications

Technical

Gross volume	296 litres
Net volume	284 litres
Temp. range	-60°C to -86°C
Temp. probes	1/PT 1000
Compressor	Embraco
Refrigerant	Nature R
Noise level	55 dB
Power consumption 24/h	8,1 kWh
Power consumption	2.957 kWh/year
Voltage	220-240V
Frequency	50 Hz

Features

Audible & visual alarm	YES
Castors	4
Integrated lock	YES
Inner lids	3
Defrost	Manual
Battery Backup	48h
USB	NO
Data connection	NO
Digital controller with display	XR30CX
Data logging	NO
High/Low temperature alarm	YES
Porthole for external probe & CO2 backup	Ø 12 mm
Prepared for external logging equipment	YES
Connection for remote alarm	YES

Material

Inner	Precoated white steel
Outer	Galvanized pre-coated steel

Outer Dimensions

Height (with castors)	885 mm
Width	1265 mm
Depth (incl. condenser)	630 mm
Depth (incl. condenser & handle)	680 mm
Net weight	76 kg

Inner dimensions

Height	635/385 mm
Width	1100/890 mm
Depth	440 mm
Insulation	80 mm
Insulation type	PUR
Internal humidity	< 5 %

ULT C300

Product no.

N113003



Storage capacity

Box height	No. of boxes	No. of racks	Total no. of boxes	Product no.
50 mm	6	3	18	N311200
	11	18	198	N311220
			= 216	
75 mm	4	3	12	N311300
	8	18	144	N311320
			= 156	
100 mm	3	3	9	N311400
	5	18	90	N311420
			= 99	

Chest freezer ULT C300

Dimensions (mm)

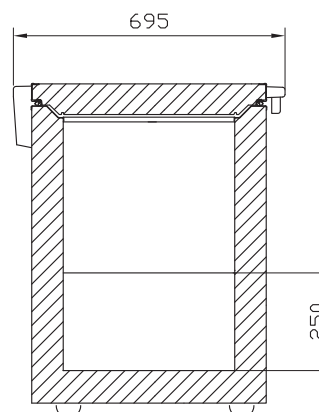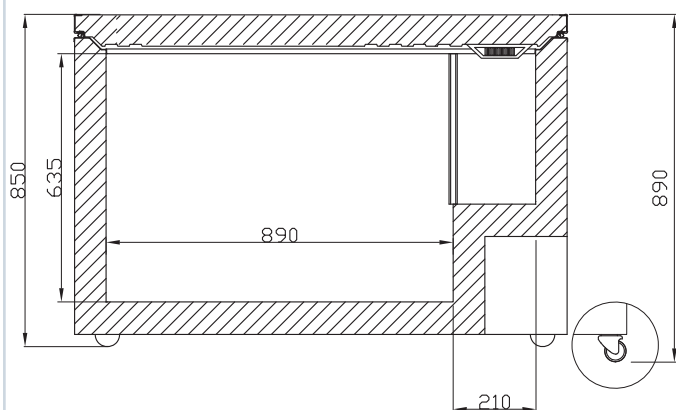ULT C300

Racks for ULT C300

ORGANIZE AND OPTIMIZE THE FREEZER

Quality, efficiency and capacity at the best possible economy, is a vital objective in Nordic Lab's business philosophy. Our storage solutions include racks and boxes, all insuring an optimal handling and market low cost per sample to be stored.

Please find below specifications on the storage solution for the ULT C300 freezer model.

Rack

ULT C300

Storage capacity

Box height	No. of boxes	No. of racks	Total no. of boxes	Racks without B&D* Product no.	Racks with B&D* Product no.
50 mm	6	3	18	N311200	N312200
	11	18	198	N311220	N312220
			=216		
75 mm	4	3	12	N311300	N312300
	8	18	144	N311320	N312320
			=156		
100 mm	3	3	9	N311400	N312400
	5	18	90	N311420	N312420
			=99		

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Warning

As the appliance contains a flammable refrigerant, it is essential to ensure that the refrigerant pipes are not damaged.

Standard EN378 specifies that the room in which you install your appliance must have a volume of 1m³ per 8 g of hydrocarbon refrigerant used in the appliances. This is to avoid the formation of flammable gas/air mixtures in the room where the appliance is located in the event of a leak in the refrigerant circuit. The quantity of the refrigerant used in your appliance is indicated on the rating plate.

WARNING:

Keep ventilation openings in the appliance's cabinet or in the built-in structure clear of obstruction

WARNING:

Do not use other **mechanical devices** or other means to accelerate the defrosting process than those recommended by the manufacturer

WARNING:

Do not damage the refrigerant system

WARNING:

Do not use **electrical appliances** inside the refrigerated storage compartment, unless they are of a type recommended by the manufacturer

WARNING:

Do not expose the appliance to rain, and secure not splashing water when cleaning the floor.

WARNING:

This appliance is not intended for use by persons (including children) with reduced physical, sensory or mental capabilities, or lack experience and knowledge, unless they have been given supervision or instructions concerning use of the appliance by a person responsible for their safety. Children should be supervised to ensure that they do not play with the appliance.

WARNING:

Children must not play with, on, or around the appliance.

WARNING:

Children must not clean the appliance or carry out general maintenance unless they are at least 8 years old and are being supervised.

WARNING:

Danger risk of fire or explosion. Flammable refrigerant used. To be repaired only by trained personnel.

**WARNING:**

Do not store explosive substances such as aerosol cans with a flammable propellant in this appliance.

WARNING:

Sharp edges on cabinet, compressor compartment, evaporator, ventilation cover and on internal equipment can occur. Please be aware to avoid injury.

WARNING:

The condenser on the back of the appliance will in some cases

have a hot surface. Please be aware to avoid injury.

- Always keep the keys in a separate place and out of reach of children
- Before servicing or cleaning the appliance, unplug the appliance from the mains or disconnect the electrical power supply
- If the supply cord is damaged, it must be replaced by the manufacturer, its service agent, or similarly qualified persons in order to avoid a hazard
- **Relevant for Australia:** Supply cord fitted with a plug complies with AS/NZS 3112.
- Frost formation on the interior evaporator wall and upper parts is a natural phenomenon. Therefore, the appliance should be defrosted during normal cleaning or maintenance.
- Please note that changes to the appliance construction will cancel all warranty and product liability

Contents

Warning2

Electrical connection.....3

Get to know your chest freezer5

Installation and start-up6

Operation and function7

Defrosting8

Cleaning and maintenance.....9

Fault finding10

Warranty, spare parts and Service11

Disposal.....12

Electrical connection

Wiring and connections in power supply systems must be all applicable (local and national) electrical codes. Consult these codes lengths and sizes prior to cabinet installation.

This device complies with relevant EU directives including Low Voltage Directive 2014/35/EU and Electromagnetic Compatibility Directive 2014/30/EU

The socket should be freely accessible.

Connect the appliance only to 220/240 V / 50Hz alternating current via a correctly installed earthed socket.

The socket must be fused with a 10 A - 13 A fuse.

If the appliance is to be operated in a non-European country, check on the rating plate whether the indicated voltage and current type correspond to the values of your mains supply.

Information regarding voltage, current or power are given on the rating plate.

The power cord may be replaced by a technician only.

The rating plate provides various technical information as well as type and serial number.

Get to know your chest freezer

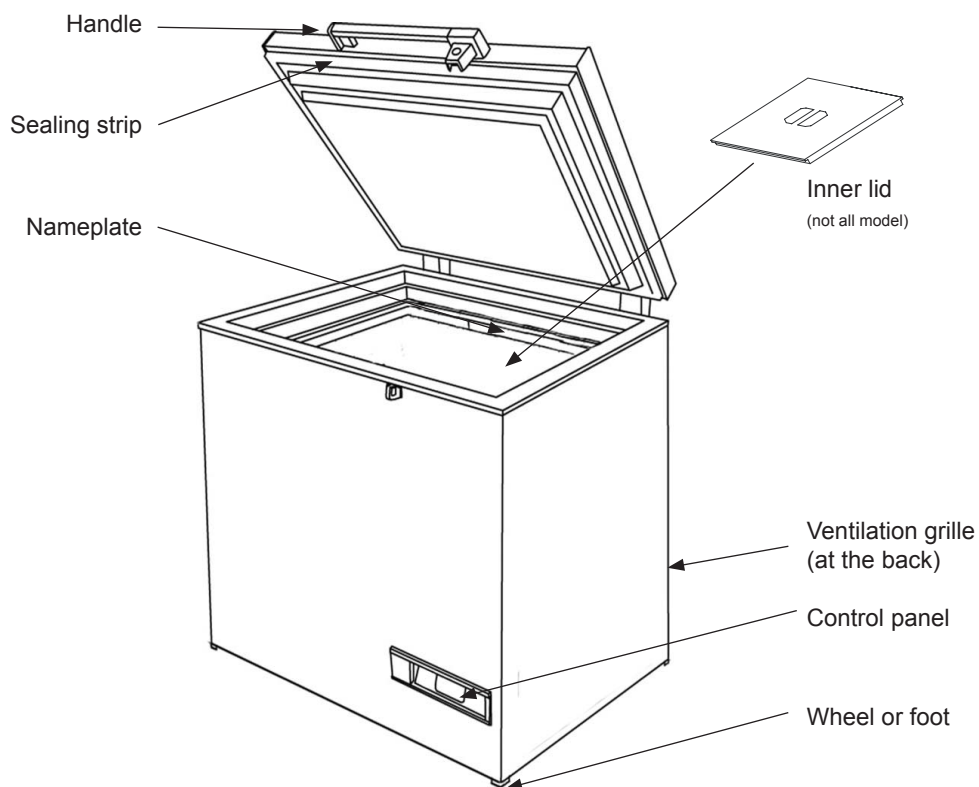
fig. 1

Installation and start-up

Room temperature

The climate class is stated on the nameplate (see page 10). This specifies the optimum room temperature.

SP = -82	SP = -86
+16 °C to +25 °C	+16 °C to +20 °C

Ventilation

It is important that the appliance be well ventilated and that air can circulate unhindered above, below and around it. Figure 2 illustrates how the necessary air circulation can be ensured.

Placement

For safety and operational reasons, the appliance must not be installed outdoors.

The appliance should be placed on a stable, level surface in a dry, well-ventilated room (max. 75% relative air humidity).

Never place the appliance close to sources of heat such as radiators, and avoid placing it in direct sunlight.

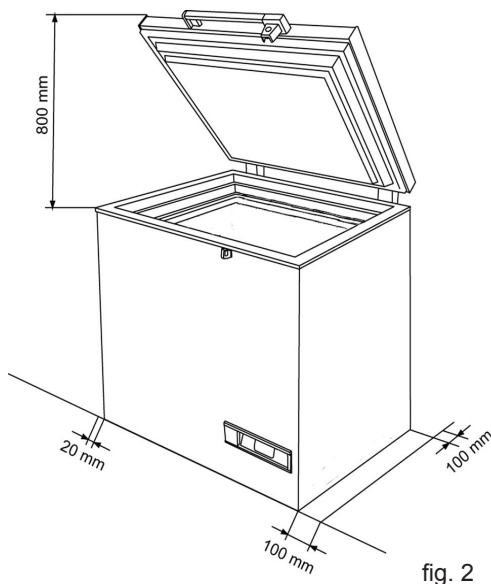
fig. 2

There must be at least 35 mm clearance between the base of the appliance and the floor.

IMPORTANT! There must be free access to the ventilation grille (fig. 3).

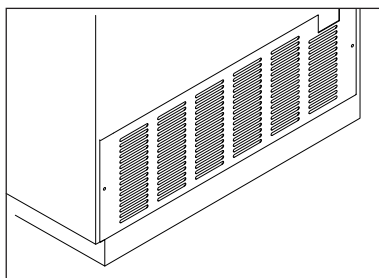
fig. 3

Operation and function

The appliance is equipped with electronic controls (fig. 4).

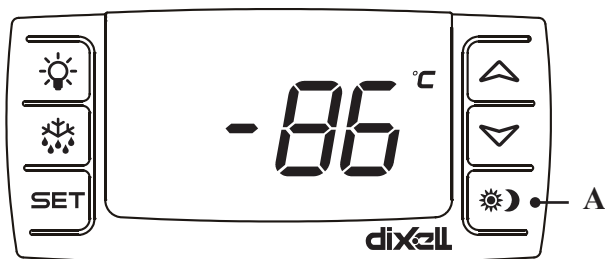
fig. 4

For operation of the electronic control, reference is made to the separate instruction manual enclosed in the appliance.

Additional information

NOTE: The first time the power is connected, press the “A” (Fig. 4), thereby the alarm will be connected


The appliance is equipped with a battery back-up. The alarm will automatically go off by power supply failure.

Therefore it is important to disconnect the alarm, in case the appliance is not to be used for a period.

Set the alarm off by pressing the “A” (fig. 4), and display shows “OFF”, and then disconnect the power supply.

Port hole

Some models has a predrilled hole at the back side. Useful in case CO₂ back up or introduction of wire sensor may be required.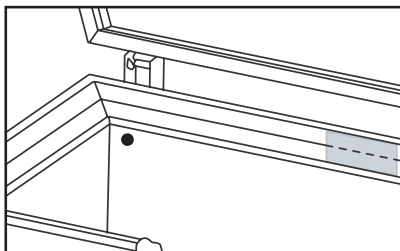

Not all models

Defrosting

The items stored in the freezer and the humidity of the surrounding air gradually cause frost and ice to form inside the freezer. A thin layer of frost or ice does not affect freezer performance, but the thicker the layer, the higher the energy consumption. Small amounts of loose frost can be removed using a plastic scraper.

Never use electrical apparatus or sharp implements as these may damage the inside of the appliance.

When the layer of frost and ice exceeds 4-5 mm the freezer should be completely defrosted. It is recommended that defrosting be performed when there are no or only a few items stored in the freezer.

Proceed as follows:

1. Pull out the power plug
2. Place a bowl of hot (but not boiling) water in the freezer, close the lid and wait approx. 30 minutes before removing loose frost with a plastic scraper.
3. When all the frost and ice have melted, clean the inside of the appliance using water with a mild, perfume-free detergent. Dry the freezer with a soft cloth.
4. Reconnect the power supply.
5. Return items to the freezer when the required temperature has been reached.

Cleaning and maintenance

Cleaning

Before cleaning, disconnect the power supply and remove all loose fittings.

The appliance is best cleaned using lukewarm water with a little mild, perfume-free detergent. Never use cleaning agents that scour. Use a soft cloth. Rinse with clean water and dry thoroughly. It is important to prevent water from entering the control panel.

The sealing strip around the lid must be cleaned regularly to prevent discolouration and prolong service life. Use clean water. After cleaning the sealing strip, check that it continues to provide a tight seal.

It is recommended that the wire and tubular condensers on the rear of the appliance be cleaned.

The ventilation grille must also be kept free of dust and dirt. Clean it with a vacuum cleaner.

If the appliance is not to be used for any length of time, switch it off, disconnect the power supply, empty it, clean the inside and leave the lid open to allow air circulation and prevent smells.

After each opening of both the lid and lower lid it is recommended to clean both the lid, the lower lid and the frame for water, condensation or the like.

Save energy

- Keep the lid closed as much as possible.
- Check regularly that the lid closes tightly.
- Position the freezer in a cool (min. 5°C), well-ventilated room.
- Defrost the freezer before the layer of ice becomes excessively thick.

Fault finding

Fault	Possible cause	Remedy
There is no light in the display	The appliance is switched off. Power failure; the fuse is blown; the appliance is not plugged in correctly.	Check the plug and socket. Check that power is connected. Reset the fuse.
Temperature in freezer too high.	The ventilation grille is blocked. The lid is not closed properly. The quantity of items placed in the freezer at a time is greater than the capacity of the appliance. Thick layer of frost and ice. The thermostat is set too high. The lid is prevented from closing tightly by items stored in the freezer.	Ensure unhindered air circulation. Close the lid and wait a few hours. Wait a few hours. Defrost the freezer. Set the thermostat to a lower temperature. Reposition items inside the freezer and check that the lid closes tightly.
Temperature in freezer too low.	The thermostat is set too low.	Set the thermostat to a higher temperature.
The freezer produces large quantities of frost.	Air humidity in the room is high. Items placed in the freezer are giving off moisture.	Ventilate the room more often or move the freezer to a room with lower humidity. Repack items stored in the freezer.
The lid cannot be opened.	Underpressure in freezer.	Wait 2-3 minutes until the underpressure has been equalised and try again.
Vibration and unusual noise.	The appliance is not on a level surface. The appliance is not free standing. Partition are loose.	Position the freezer on a flat, level and stable surface. Move the appliance away from other appliances or pieces of furniture. Check movable fittings and reposition them.

Warranty, spare parts and service

Warranty disclaimer

Faults and damage caused directly or indirectly by incorrect operation, misuse, insufficient maintenance, incorrect building, installation or mains connection. Fire, accident, lightening, voltage variation or other electrical interference, including defective fuses or faults in mains installations are not covered by the warranty.

Repairs performed by others than approved service centres and any other faults and damage that the manufacturer can substantiate are caused by reasons other than manufacturing or material faults are not covered by the warranty.


Please note that changes to the construction of the appliance or changes to the component equipment of the appliance will invalidate warranty and product liability, and the appliance cannot be used lawfully. The approval stated on rating plate will also be invalidated.

Transport damage discovered by the buyer is primarily a matter to be settled between the buyer and the distributor, i.e. the distributor must ensure that such complaints are resolved to the buyer's satisfaction.

Before calling for technical assistance, please check whether you are able to rectify the fault yourself. If your request for assistance is unwarranted, e.g. if the appliance has failed as a result of a blown fuse or incorrect operation, you will be charged the costs incurred by your call for technical assistance.

Spare parts

When ordering spare parts, please state the type, serial and product numbers of your appliance. This information is given on the rating plate. The rating plate contains various technical information, including type and serial numbers.

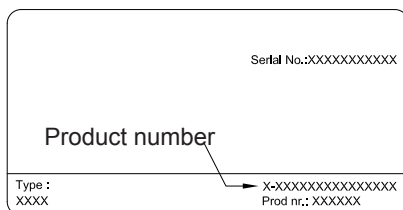Disposal

Information for Users on Collection and Disposal of Old Equipment and used Batteries

These symbols on the products, packaging, and/or accompanying documents mean that used electrical and electronic products and batteries should not be mixed with general household waste. For proper treatment, recovery and recycling of old products and used batteries, please take them to applicable collection points, in accordance with your national legislation and the Directives 2012/19/EU and 2006/66/EC.

By disposing of these products and batteries correctly, you will help to save valuable resources and prevent any potential negative effects on human health and the environment which could otherwise arise from inappropriate waste handling.

For more information about collection and recycling of old products and batteries, please contact your local municipality, your waste disposal service or the point of sale where you purchased the items.

Penalties may be applicable for incorrect disposal of this waste, in accordance with national legislation.

For business users in the European Union.

If you wish to discard electrical and electronic equipment, please contact your dealer or supplier for further information.

Information on Disposal in other Countries outside the European Union

These symbols are only valid in the European Union. If you wish to discard this product, please contact your local authorities or dealer and ask for the correct method of disposal.

Note for the battery symbol

This symbol might be used in combination with a chemical symbol. In this case it complies with the requirement set by the Directive for the chemical involved.

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

Maintenance for Chest Freezers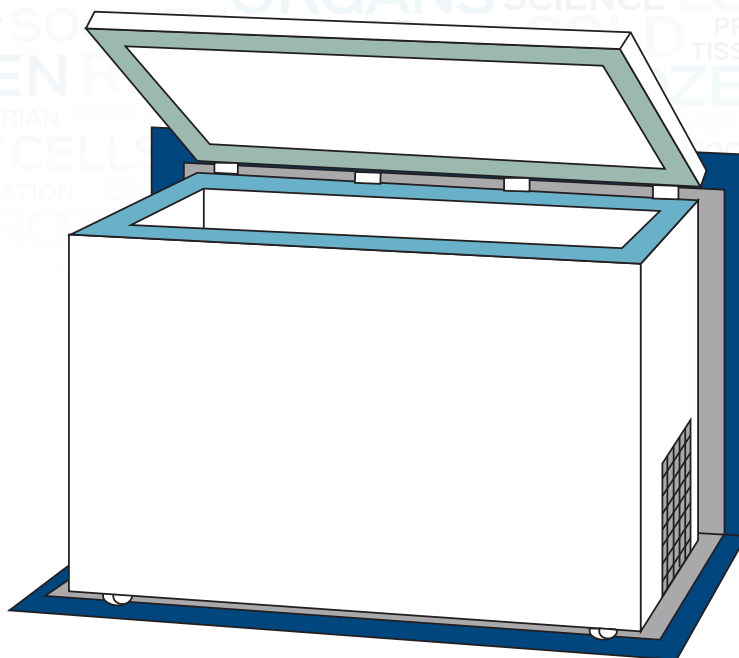

Maintenance										
Dark blue	Weekly	Check placement: Min.10 cm free space to the side of the freezer Min. 10 cm free space to the back of the freezer								
Light blue	Weekly	Defrosting: Small amounts of loose ice can be removed using a plastic scraper. Defrosting has to be done regularly to prevent ice building The ice build depends on the use of the freezer. The more daily lid openings - the more manual defrosting. <table><tr><td>Lid open/close per 24h</td><td>Advised monthly manual defrosting</td></tr><tr><td>< 5</td><td>1-2</td></tr><tr><td>< 10</td><td>2-4</td></tr><tr><td>< 15</td><td>4-6</td></tr></table>	Lid open/close per 24h	Advised monthly manual defrosting	< 5	1-2	< 10	2-4	< 15	4-6
Lid open/close per 24h	Advised monthly manual defrosting									
< 5	1-2									
< 10	2-4									
< 15	4-6									
Gray	Monthly	Clean: The ventilation grille, the condenser, the compressor, the compressor compartment and the floor must also be kept free of dust and dirt. Clean it with a vacuum cleaner regularly.								
Green	Quarterly	Clean: The sealing strip around the freezer lid must be cleaned regularly to prevent discolouration and prolong service life. Use clean water. After cleaning the sealing strip, check that it continues to provide a tight seal.								

Maintenance chart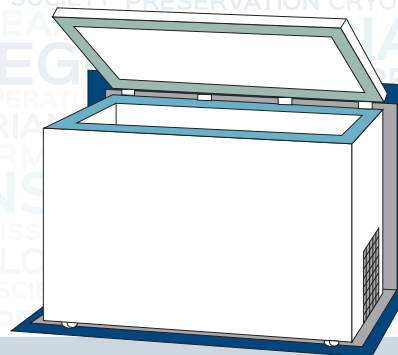

Maintenance chart

Dark blue

Weekly

Check placement:

Date:

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

Light blue

Weekly

Door openings

per 24h / Monthly

manual defrosting

< 5 / 1-2

< 10 / 2-4

< 15 / 4-6

Date:

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

Gray

Monthly

Clean:

Date:

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

Green

Quarterly

Clean:

Date:

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—