AUTOMATIC BURRETE ASSEMBLY STEP BY STEP



BURETTE WITH DETACHABLE PTFE KEY ASSEMBLY INSTRUCTIONS:


AUTOMATIC BURRETE ASSEMBLY STEP BY STEP



Assemble clamp on bottle neck, & place springs on possition



Adapter consist of center part/connector, piece of silicon sleeve, two outer caps/screw thread.

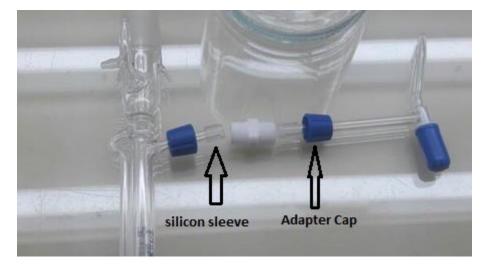


AUTOMATIC BURRETE ASSEMBLY STEP BY STEP

Insert screw cap of adapter on both ends of tubes with threads on outer side, Put silicon sleeve on both ends of inclined tube of automatic burette.



Joint both tubes with center part of adapter. Tighten screw to avoid leakage.

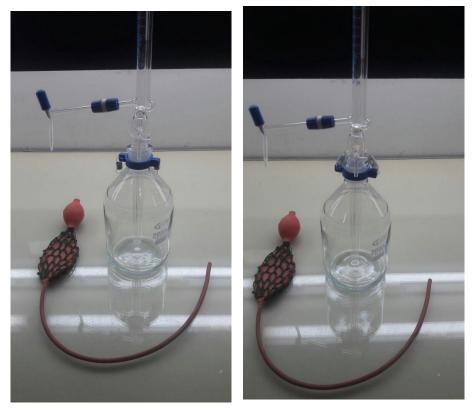

Mount burette on bottle & connect rubber ballow & springs to burette hooks.