TR06 Tube roller

Please read the User Manual carefully before use, and follow all operating and safety instructions!

user manual

english / español / français / italiano

User Manual

TR06 Tube roller, variable speed

Preface

Users should read this Manual carefully, follow the instructions and procedures, and beware of all the cautions when using this instrument.

Service

In order to guarantee this equipment Works safely and efficiently, it must have a regular maintenance. In case of any faults, do not try to repair it yourself. If help is needed, you can always contact your dealer or Labbox via www.labbox.com

Please, provide the customer service representative with the following information:

- Serial number
- Description of the problem
- Your contact information

Warranty

This instrument is guaranteed to be free from defects in materials and workmanship under normal use and service, for a period of 24 months from the date of invoice. The warranty is extended only to the original purchaser. It shall not apply to any product or parts which have been damaged on account of improper installation, improper connections, misuse, accident or abnormal conditions of operation.

For claim under the warranty, please contact your supplier.

Warning: do not use the rotor if it has been damaged due to corrosion and/or fatigue.

1. Safety Instructions

	<p>Warning!</p> <ul style="list-style-type: none">• Read the operating instructions carefully before use.• Ensure that only trained staff works with the instrument.
	<p>Protective ground contact!</p> <ul style="list-style-type: none">• Make sure that socket must be grounded (protective ground contact) before use.

Table 1

- When working, wear the personal guard to avoid the risk from:
 - Splashing liquids
 - Broken glass containers
- Follow the safety instructions, guidelines and accident prevention regulations.
- Do not touch the running parts, moving instrument care not rolling your fingers.
- Set up the instrument in a spacious area on a stable, clean, non-slip, dry and fireproof surface. Do not operate the instrument in explosive atmospheres, with hazardous substances or under water.
- If the instrument does not run smoothly, please decrease the motor speed.
- Preparation of samples may lead to dangerous flammable. Only process samples that will not react dangerous.
- The instrument only be opened by expert, please switch off before use.
- The voltage stated on the nameplate must correspond to the mains voltage.
- Do not cover the instrument during running. Prevent the collision and extrusion to instrument and accessories.

2. Proper Use

The instrument is designed for mixing substances in schools, laboratories or factories. This device is not suitable for using in residential areas or other constraints mentioned in Chapter 1. Failure to use the instructions may be caused unsafe situation.

3. Inspection

3.1. Receiving Inspection

Unpack the equipment carefully and check for any damages which may have arisen during transport. If it happens, please contact manufacturer/supplier for technical support.

Note:
If there is any apparent damage to the system, please do not plug it into the power line.

3.2. Listing of Items

The package includes the following items:

Items	Qty
Main unit	1
Power cable	1
User Manual	1

Table 2

4. Control

Figure 1

Items	Default settings
Speed control knob	The rotating function is switched ON/OFF by rotating the knob. To increase the value turn the knob clockwise.
Power light Power	Green LED display light. LED light shows green when switch on.
Power switch I/O	Switch ON or OFF the instrument.

Table 3

5. Operation

- Place the Rotator in safe and stable surface, ensure the required operating voltage and power supply voltage matched, and the socket must be earthed reliably.
- Connect the power cable.
- Switch on instrument by pressing the power switch I/O.
- Rotate clockwise the speed control knob to start rotating function and to set target speed.
- At the end of the manipulation, turn the knob fully counterclockwise to turn off the rotation.
- Turn off the instrument by pressing the power switch I/O.

If these operations above are normal, the instrument is ready to operate. If not, the instrument may be damaged during transportation, please contact technical support of manufacturer/supplier.

Note:

It is forbidden to manipulate / move the container when the unit is in operation.
The set values can be altered at any time.

6. Faults

- Instrument cannot be powered ON.
 - Check whether the power cable is connected.
- Rotating function suddenly stop
 - Turn the knob fully counterclockwise to turn off the rotation, and then restart rotating function after lowered load. If rotating function cannot start, then repeat the process and gradually reduce load.

If these faults are not resolved, please contact manufacturer/supplier.

7. Maintenance and Cleaning

- Proper maintenance can keep instruments working properly and lengthen its lifetime.
- Do not spray cleanser into the instrument when cleaning.
- Remove the power line when cleaning.
- Only use recommended cleansers:

Dyes	Isopropyl alcohol
Construction materials	Water containing tenside / Isopropyl alcohol
Cosmetics	Water containing tenside / Isopropyl alcohol
Foodstuffs	Water containing tenside
Fuels	Water containing tenside

Table 4

- Before using other method for cleaning or decontamination, the user must ascertain with the manufacturer that this method will not harm the instrument.
- Wear the proper protective gloves during cleaning of the instrument.

**Note:**

- Electronic device cannot clean with cleanser.
- If you require maintenance service, must be cleaned the instrument in advance to avoid pollution of hazardous substances, and to send back into original packing.
- If the instrument will not use for a long time, please switch off and place in a dry, clean, room temperature and stable location.

8. Associated Standards and Regulations

Construction in accordance with the following safety standards:

EN 61010-1

UL 3101-1

CAN/CSA C22.2(1010-1)

EN 61010-2-10

Construction in accordance with the following EMC standards:

EN 61326-1

Changes or modifications not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment.

9. Specifications

Items	Specifications
	TR06 series
Voltage [VAC]	100-240
Frequency [Hz]	50/60
Power [W]	25
Number of rollers	6
Roller size [length x diameter, mm]	280 x 30
Motor type	DC motor
Speed range [rpm]	0-70
Run type	Continuous operation
Overall Dimensions [mm]	470×260×120
Weight [kg]	4.5
Permissible ambient temperature [°C]	5-40
Permissible relative humidity	80%
Protection class acc. to DIN EN60529	IP21

Table 5

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