## **ULTR**      **Ultrasonic bath with heater and timer**

*Please read the User Manual carefully before use, and follow all operating and safety instructions!*

# **user manual**

english / español / français / italiano

## User Manual

# Ultrasonic bath with heater and timer

## Preface

Users should read this Manual carefully, follow the instructions and procedures, and beware of all the cautions when using this instrument.

## Service

When help is needed, you can always contact the Service Department of manufacturer for technical support:

**[www.labbox.com](http://www.labbox.com) / e-mail: [info@labbox.com](mailto:info@labbox.com)**

Please provide the customer care representative with the following information:

- Serial number
- Description of problem
- Your contact information

## Warranty

This instrument is warranted to be free from defects in materials and workmanship under normal use and service, for a period of 24 months from the date of invoice. The warranty is extended only to the original purchaser. It shall not apply to any product or parts which have been damaged on account of improper installation, improper connections, misuse, accident or abnormal conditions of operation.

For claim under the warranty please contact your supplier.

# 1. Safety Instructions

	<b>Warning!</b> <ul style="list-style-type: none"><li>• Read the operating instructions carefully before use.</li><li>• Ensure that only trained staff works with the instrument.</li></ul>
	<ul style="list-style-type: none"><li>• <b>Do not operate if the tank is empty</b>, or the ultrasonic generator will be damaged. If heating is needed, <b>the water level shouldn't be less than 2/3.</b></li></ul>
	<ul style="list-style-type: none"><li>• <b>Do not turn on the machine if the tank is empty or the transducer will be damaged</b></li><li>• Do not relocate the machine when there is fluid in the tank (to avoid overflow)</li><li>• Water-soluble liquids are suggested for the ultrasonic cleaner. <b>Strong acids or flammable cleaners are strictly forbidden.</b> If a solvent type cleaner is added in the tank, please stop working for a while after the 30-minute cycle</li></ul>
	<ul style="list-style-type: none"><li>• Connect the device to an earthed power supply to ensure the safety of the machine and the experiment; connect the power as the machine requires it.</li><li>• Refitting is strictly prohibited!</li><li>• Ensure the earth wire is well connected before starting.</li></ul>

Table 1

- The control panel cannot be cleaned with organic solutions, strong acids or strong alkalis it could become eroded or damaged.
- Please close the lid to reduce noise and pay attention to the water and the steam they could cause skin burns.

### **Do not use the machine in any of the environments below:**

- A place where temperature changes fiercely
- A place where vibration or impact is strong
- A place where water oil or chemicals splash
- A place where humidity is too high and it is easy to produce dew
- A place where there is corrosive gas or dust
- A place where there is explosive and flammable gas

## 2. Inspection

### 2.1 Receiving Inspection


Unpack the equipment carefully and check for any damages which may have arisen during transport. Please contact your supplier for technical support.

**Note:**

If there is any apparent damage to the system, please do not plug it into the power line.


### 2.2 Specifications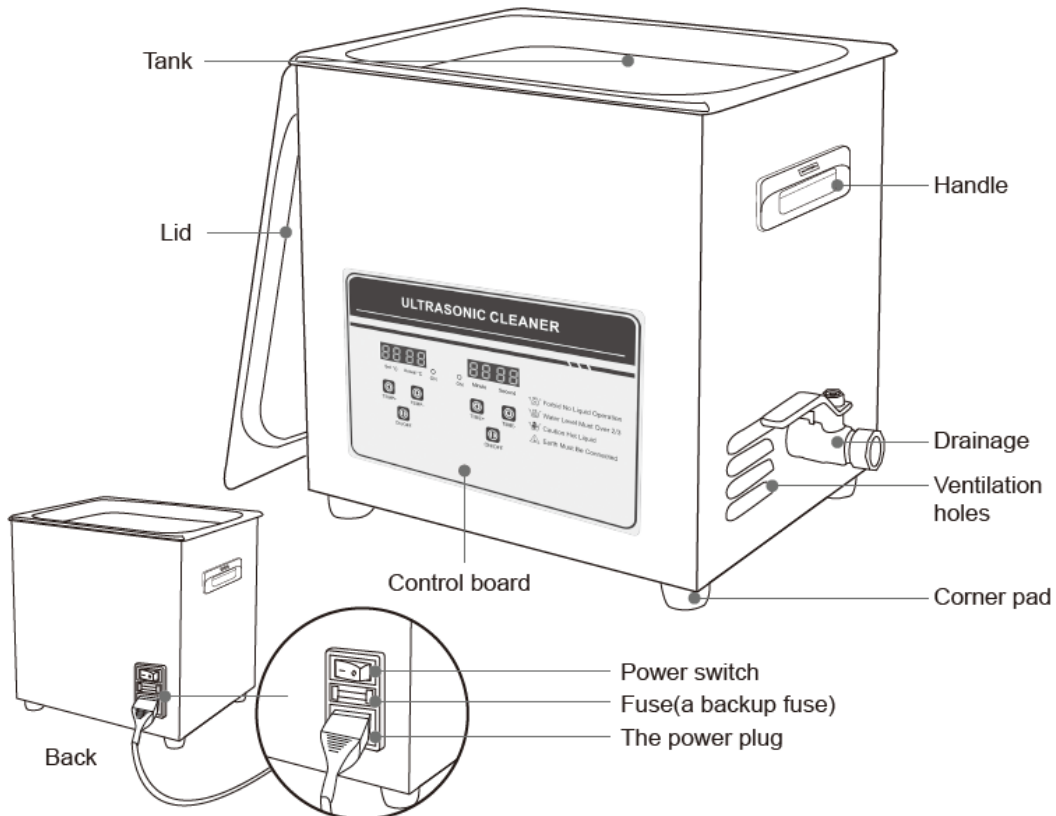

**\*Drainage only available in 6,5 L – 30 L models**


Model	Frequency	Transducer (pcs)	Capacity (L)	Tank size (mm)	Unit size (mm)	Heating Power (W)	Ultrasonic Power (W)	Timer (M)	Temp. (°C)
ULTR-2L0-001	40	1	2	150x135x100	175x160x210	100	80	0-30	RT-80
ULTR-3L2-001	40	2	3.2	240x135x100	265x165x220	100	120	0-30	RT-80
ULTR-6L5-001	40	3	6.5	300x150x150	325x180x280	200	180	0-30	RT-80
ULTR-10L-001	40	4	10	300x240x150	325x265x280	200	240	0-30	RT-80
ULTR-15L-001	40	6	15	330x300x150	360x325x285	300	360	0-30	RT-80
ULTR-30L-001	40	10	30	500x300x200	530x325x325	500	600	0-30	RT-80

### 3. Operation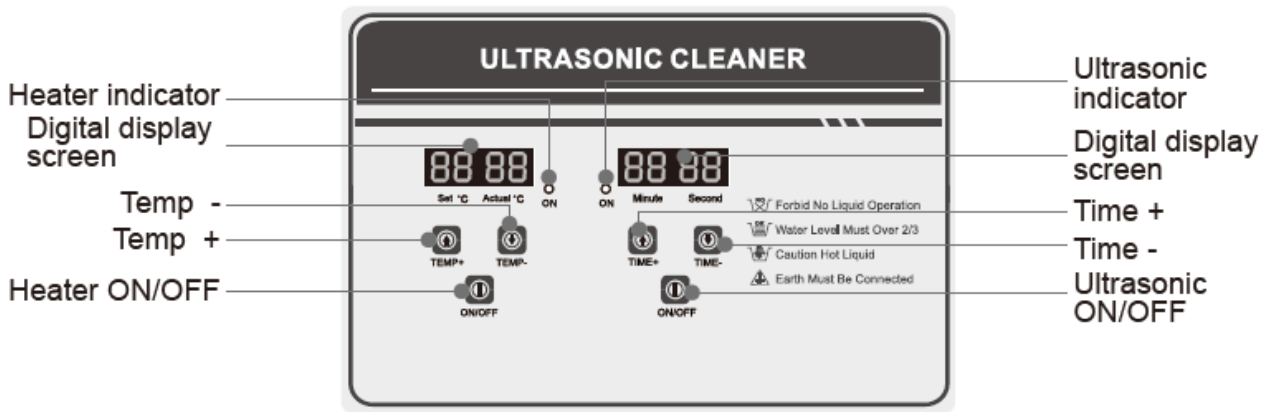


#### Operation steps:

##### ◆ Before operating your cleaner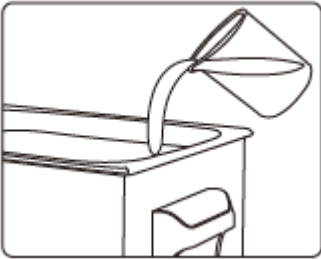

##### 1. Add cleaning liquid

Fill the tank with warm or tap water below the maximum line; Add the required quantity of cleaning solution. When using flammable liquid such as alcohol, gasoline and so on, heating is strictly forbidden! Ultrasonic vibration for a long time can cause temperature of liquid increase, caution fire!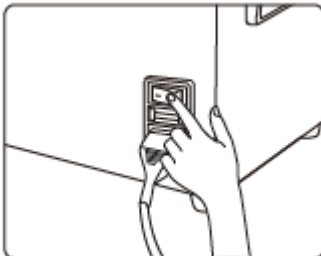

##### 2. Turn on the power switch


Plug the cleaner into a grounded outlet. Turn on the power switch on the back of the tank. ("O" is closed, and "I" is open)

##### ◆ Start ultrasonic cleaning

For the first boot, time display show setting time at random, press **time key+** one time, "the setting time" on the display will be increase by 1 minute correspondingly, touch **time key-** to decrease time setting by 1 minute. Touch **ON/OFF** switch after you have properly set the time, ultrasonic indicator will turn red and the machine will start working based on the parameters set. Meanwhile the timer will begin to count down and stop working when the time shows "00:00", automatically.

## ◆ Heat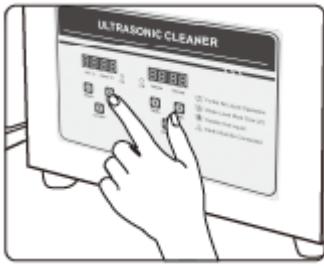

For the first boot, the setting temperature on the display is random, and actual temperature is room temperature. Touch **TEMP+** once to increase temperature by 1°C; Touch **TEMP-** once to decrease temperature by 1°C. Temperature display shows setting temperature and actual temperature when heater is working, the heater indicator goes off when actual temperature reaches setting temperature.

## ◆ Power OFF

Press “ON/OFF” to stop working and then cut off the power. Pour the liquid out, clean the tank and the unit with cloth for next use.

## 4. Maintenance

- To keep your device in good working conditions the electrical control system should be blown regularly with high pressure air to get rid of the dust.
- Clean the tank at least once a week

### Trouble clearing

No.	Problems	Possible Cause	Solutions
1	No ultrasonic activity	A. Power supply not connected B. Fuse broken C. Cable is short circuited D. Transducer is short circuited E. PCB Board broken F. Other reasons	A. Check and plug power switch B. Check the fitted power supply and change the same specification fuse C. Connect fitted cable or replace a new D. Consult with your supplier E. Check the broken part and change it F. Consult with your supplier
2	Time control failure	A. Timer failure B. Other reasons	A. Replace the timer or digital control panel B. Consult with your supplies
3	No heating	A. Heater power switch had a bad linkage B. Fuse burn out C. Heating pad burn out D. Digital display board out of control E. Other reasons	A. Check heating plug and connect it well. Check outlet line with multimeter: OK if resistance value is a few hundred OHMs. B. Change the same specification fuse C. Replace the bad heating pad if open circuit D. Check the broken part and change it E. Consult with your supplier
4	Temperature control failure	A. Thermostat loosen B. Thermostat tube broken C. Digital display out of control D. Other reasons	A. Fasten the thermostat header B. Replace the thermostat C. Check the broken part and change it D. Consult with your supplier
5	Not well cleaning	A. Not strong ultrasonic cleaning B. Too high or too low liquid surface C. Too high too low temperature D. Not suitable cleaning liquid E. Other Reasons	A. Connect ultrasonic button and adjust B. Adjust liquid into the best surface C. Adjust the temperature into the most fitted D. Stop and switch off power supply, replace suitable liquid after previous liquid cooling down E. Consult with your supplier
6	Electric Leakage	A. Customer side not grounded B. Machine not grounded	A. Ensure it is grounded B. Check if machine earth wire loosens
7	Scrap		Not repairable or discarded

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