





## Product datasheet

### Captair 391 Smart


Secure weighing station

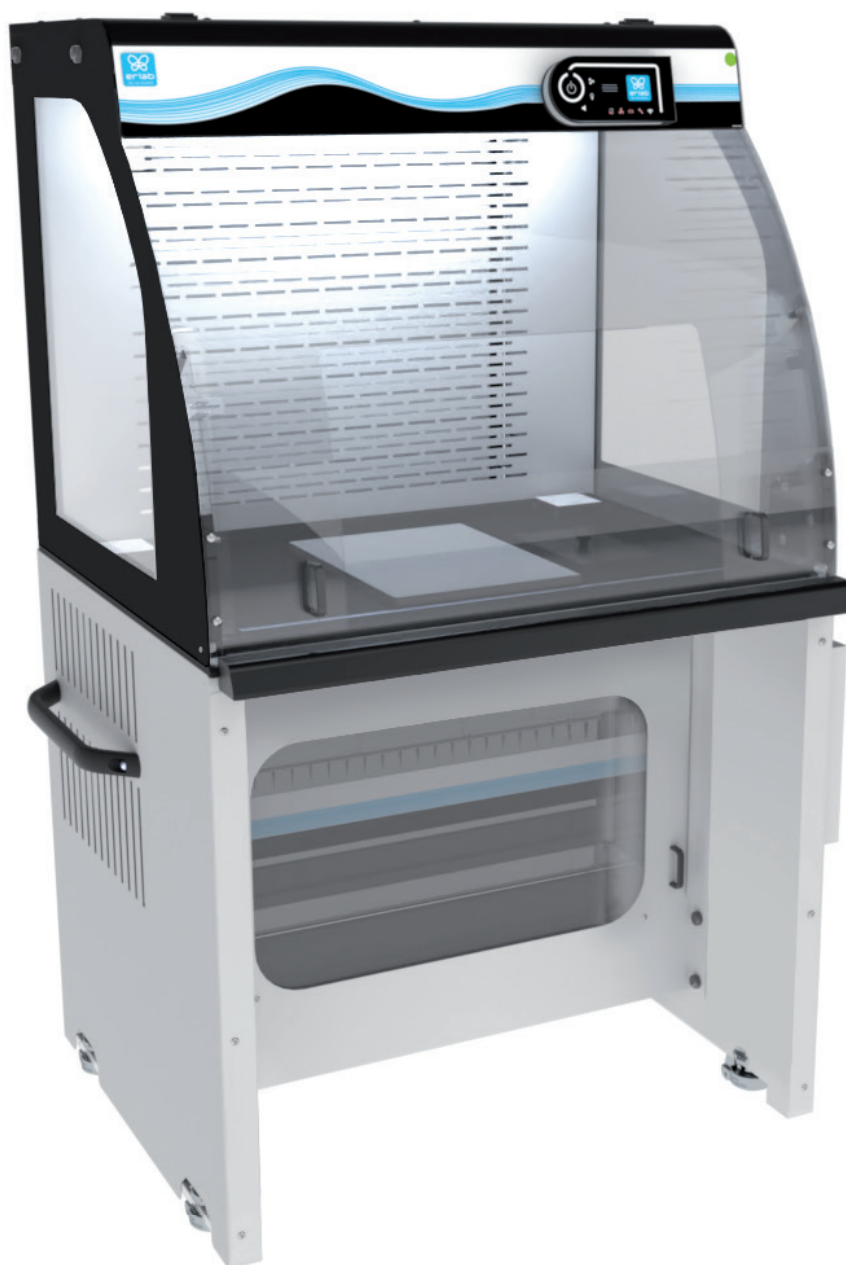

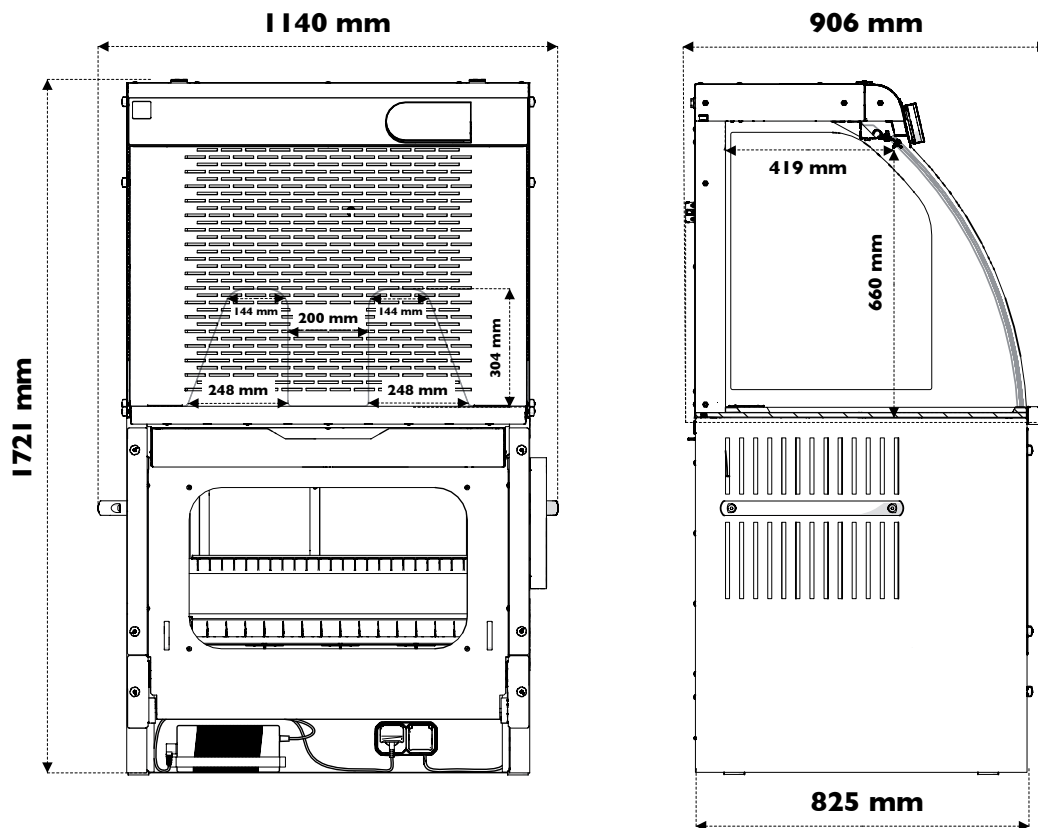

Work surface with built-in spill tray

**Trespa® Top Lab<sup>PLUS</sup>**

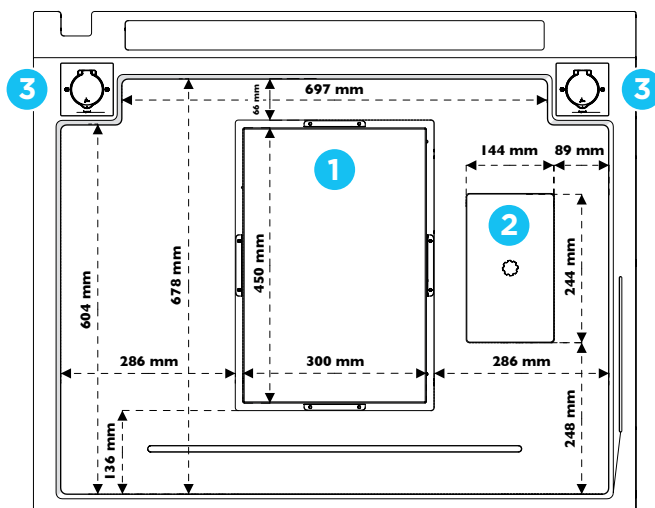




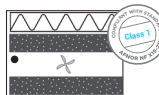
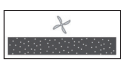
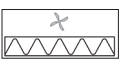
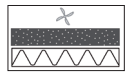
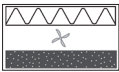
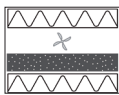
- ① Weighing plate
- ② Waste port
- ③ Electrical outlet

Fixed work bench

## FILTRATION TECHNOLOGY



Modular design of the filtration column allows to adapt to every protection needs.

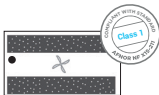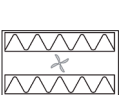		Products handled / Applications			
		Liquid chemicals handlings	Powders handlings	Liquid chemicals and powders handlings	Liquid chemicals handlings in clean room
Customized filtration column	Class 1 according to the NF X15-211 standard	 <b>2C</b>	 <b>2P</b>	 <b>1P2C</b>	 <b>2C1P</b>
	Class 2 according to the NF X15-211 standard	 <b>1C</b>	 <b>1P</b>	 <b>1P1C</b>	 <b>1C1P</b>  <b>1P1C1P</b>



### Carbon filtration for gases and vapours

AS: For Organic vapours  
BE+: Polyvalent for Acid + Organic vapours  
F: For Formaldehyde vapours  
K: For Ammonia vapours



### Particulate filtration for powders

HEPA H14: 99.995% efficiency filtration of particles over 0.1µm in size  
ULPA U17: 99.99995% efficiency filtration of particles over 0.1µm in size



- **Molecode**  
Automatic alarm to detect a filtration fault



**Class 1**  
= **Maximum safety**

<b>Safety standards</b>	AFNOR NF X15-211: 2009: France – BS 7989: England DIN 12 927: Germany – EN 1822: 1998 (HEPA H14 Filter) – EU Marking
<b>Air flow</b>	220m³/h (Carbon Filter) – 300m³/h (HEPA Filter)
<b>Air face velocity</b>	0.4 to 0.6m/s
<b>Voltage/Frequency</b>	110-230V/50-60Hz
<b>Power consumption</b>	Max. 2300W (with 2 sockets inside)
<b>Sash opening</b>	Oblong (Carbon Filter) or Trapezoidal (HEPA Filter)
<b>Structure</b>	Corrosion resistant electro-galvanized steel coated with antiacid polymer
<b>Side and front panels</b>	Chemical resistant acrylic
<b>Filtration module</b>	Polypropylene

## Features

<b>Communication interface</b>	Simple communication by audible and light pulses: unit running time, air face velocity, automatic alarm to detect a filtration fault, ventilation settings, fan failure alarm
<b>Filtration technology</b>	1 adaptable filtration column (with BIBO* secure filtration unit)
<b>Carbon filtration for gases and vapours</b>	Following filtration column configuration (see table above)
<b>Particulate filtration for powders</b>	Following filtration column configuration (see table above)
<b>Monitoring</b>	Real-time control of security settings
<b>Monitoring of ambient handling conditions</b>	Temperature (T°) / Hygrometry (RH) sensors
<b>Internal lighting</b>	LED lighting > 650lux
<b>Anemometer</b>	Air face velocity alarm / Air face velocity indicator
<b>Chemical Listing</b>	List of 700+ approved chemicals compliant with AFNOR NF X15-211 filtration standards
<b>Ceiling lighting</b>	ON/OFF light button
<b>Work surface</b>	Trespa® Top Lab <sup>PLUS</sup>
<b>Bench</b>	Mobile (installation) and Fixed (with anti-vibration rubber-tyred wheels)

## Options

<b>Molecode</b>	Detection sensor: Type A, for Acids / Type F, for Formaldehyde / Type S, for Solvents
-----------------	---

\*BIBO: Bag In Bag Out system.



# About ERLAB

*The ERLAB Research and Development Laboratory*

Since 1968, **ERLAB** has been a specialist, inventor and world leader in **ductless, zero-emission filtering fume hoods for laboratories** to provide total safety in chemical handling.

## 1 ERLAB filtration

We provide technologies to protect laboratory staff from inhaling chemicals. This is made possible thanks to our **Research and Development (R&D) department**, which has continuously improved our filtration technology **for more than 50 years**. That's why, in 2009, we invented the **ERLAB ABOVE** label for tried and tested filtration technology.

## 2 The AFNOR NF X15-211: 2009 standard

ERLAB's filtration technology conforms to the **NF X15-211: 2009 standard**, the industry's most demanding standard for molecular filtration, developed by a committee of independent scientists and specialized manufacturers.

**This text imposes performance criteria linked to:**

- Filtration efficiency
- Containment efficiency
- Air face velocity
- Documentation: **chemical listing**

## 3 The ESP programme

A set of three services included with the purchase of each device designed to ensure your safety.



**eValiQuest** Risk analysis – Determination of protection needs – Determination of ergonomic needs



**ValiPass** Certified installation – Total safety for handling



**ValiGuard** Ongoing monitoring – Preventative and maintenance inspections – Device reconfiguration based on protection needs – Development of handling

## 4 Flex technology

The combination of molecular and particulate filtration technologies allows a single device to meet laboratories' protection needs. This innovation from ERLAB's R&D department offers unprecedented **flexibility, versatility and value**. A single device can be reconfigured over time and easily reassigned to other applications.

## 5 Smart technology

Smart technology is a **simple and innovative** means of communication that improves safety. This technology uses a light and sound signal to indicate the user's level of protection. The advantages of the technology are:

- 1 | **Light pulsation:** Real-time communication via **LED light pulses** intuitively alerts the user to the device's operating status.
- 2 | **Simplicity:** One-touch activation.
- 3 | **Detection system:** The exclusive detection system continuously monitors filtration performance.
- 4 | **Built-in monitoring:** This service provides direct access to the **status, settings and history** of your device.

BL

**BLANC - LABO S.A.**

Chemin des Mouettes 1, Centre du Bief, CH - 1027 Lonay

Direct +41 (0)21 804 18 56

Fax +41 (0)21 804 18 59 [info@blanc-labo.com](mailto:info@blanc-labo.com)

[www.blanc-labo.com](http://www.blanc-labo.com)