Operating manual

ULT | ULTEco

ULT freezer

Table of Contents

1. About this Manual	4
2. Safety	6
2.1 Terms and Symbols Used	6
2.1.1 Terms Used	6
2.1.2 Symbols Used	6
2.2 Product Safety and Dangers	6
2.3 Requirements to be met by Operating Personnel	7
2.4 Responsibility of the Owner	7
2.5 Product Use	8
2.5.1 Intended Use	8
2.5.2 Improper Use	8
2.6 Changes and Alterations	9
2.7 Behaviour in case of Malfunctions and Irregularities	9
2.8 Switching off the Unit in an Emergency	9
3. Construction and Description	10
3.1 Design	10
3.2 Description of Function	10
3.3 Materials	10
3.4 Connections and Interfaces	11
3.4.1 Electrical Connection	11
3.4.2 Communication Interfaces	11
3.4.3 Potential-free contact	13
3.5 Nameplate	14
3.6 Technical Data	14
3.7 Applied Directives and Standards	16
3.7.1 Material Compliance	16
3.8 Ambient Conditions	16
3.9 Scope of Delivery	17
4. Delivery, Transport and Setting Up	18
4.1 Safety	18
4.2 Delivery	18
4.3 Transport	18
4.4 Unpacking	18
4.5 Storage after Delivery	19
4.6 Setting Up	19
4.6.1 Preconditions	19
4.6.2 Spacer	20
5. Putting into Operation	21

5.1	Putting into Operation for the First Time	21
5.2	Connecting the Unit to the Power Supply	21
5.3	Switching on Unit	22
6.	Operation and Control	23
6.1	Operating Personnel	23
6.2	Opening the Door	23
6.3	Switching on Unit	23
6.4	Role-based access control system	23
6.5	Log in	24
6.6	Dashboard	24
6.6.1	Distance View	25
6.6.2	Adjusting temperature	25
6.6.3	Display temperature graph	25
6.7	Menu	26
6.7.1	Language settings	26
6.7.2	Switch off appliance	26
6.7.3	Transfer log data	27
6.7.4	View battery information	27
6.8	Options	28
6.8.1	Record service activities	28
6.8.2	Display temperature and compressor status	28
6.8.3	Setting Parameters	29
6.9	Defrost the unit	31
7.	Malfunctions, Warning and Error Messages	32
7.1	Fault Messages	32
7.1.1	Dealing with fault reports	32
7.1.2	Fault Messages	33
7.1.3	Door open notification	33
7.1.4	Emergency mode	33
7.1.5	Connection error	34
7.1.6	Power Failure	34
8.	Maintenance and Servicing	35
8.1	Cleaning	35
8.2	Regular Maintenance	36
8.3	Repairs and Service	36
9.	Storage, Transport and Disposal	37
9.1	Storage and Transport	37
9.2	Disposal	37

1. About this Manual

Purpose and target audience

This manual describes the design, function, transport, operation and maintenance of the product series ULT freezer. It is intended for use by trained personnel employed by the owner who are tasked with operating and/or maintaining the unit.

If you have been tasked with working on the unit, read this manual carefully before starting work. Familiarise yourself with the safety instructions. Only perform work that is described in this manual. If there is anything you do not understand, or if any information is lacking, ask your line manager or contact the manufacturer. Do not take any course of action on your own initiative.

Versions

The appliances are available in different equipment versions and sizes. If certain features or functions are only available in certain equipment versions, this is indicated at the relevant points in this manual.

The functions described in this manual relate to the most recent firmware version.

Due to the different equipment versions and sizes, the illustrations in this manual may be slightly different to your product. However, the product is identical in terms of its operation and function.

Further applicable documents

In addition to this manual, please also refer to the following documents:

- Service manual:

To carry out service and repair work you will require the separate service manual. Manuals can be requested from Memmert International After Sales or downloaded from **www.memmert.com**.

Retaining and passing on this manual

This operating manual belongs to the unit and must always be kept in a location where it can be easily found by those working with the unit. It is the responsibility of the owner to ensure that persons who work on the unit know where this operating manual is. We recommend always storing it in a safe place near the unit.

Ensure that the manual is not damaged by heat or humidity. If the unit is sold or transported and re-installed at another location, this operating manual must be handed over with the unit. The current version of this operating manual is also available in PDF format at **www.memmert.com**.

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E-mail: service@memmert.com

www.memmert.com

If you have any queries, please always quote the product number on the nameplate.

Shipping address for repairs

Memmert GmbH + Co. KG

Willi-Memmert-Straße 90-96 | D-91186 Büchenbach | Germany

Please contact our customer service before sending appliances for repair or before making returns, otherwise, we have to refuse acceptance of the shipment.

2. Safety

2.1 Terms and Symbols Used

In this manual and on the unit itself, certain recurring terms and symbols are used to warn you of hazards or give you information that is important in order to prevent injury or damage. To avoid accidents and damage, observe and follow these instructions. These terms and symbols are explained below.

2.1.1 Terms Used

	Warns of a dangerous situation that will result directly in death or serious (irreversible) injury.
	Warns of a dangerous situation that could result in death or serious physical injury.
	Warns of a dangerous situation that could result in moderate or minor physical injury.
	Warns of damage to property.

2.1.2 Symbols Used

	Gases / vapours		Risk of explosion
	Do not lift unit		Do not tilt
	Do not enter		General warning sign
	Danger of electrocution		Tipping hazard
	Disconnect the mains plug		Wear gloves
	Wear safety shoes		Observe information in separate manual

2.2 Product Safety and Dangers

The units described in this manual are technically sophisticated, manufactured using high-quality materials and subject to many hours of testing in the factory. They reflect the state of the art and comply with recognised technical safety regulations. However, there are still risks involved, even when the units are used as intended. These are described below.

⚠ DANGER**Danger of suffocation inside the appliance**

If the appliances is a certain size, you can get accidentally locked in, which is potentially life-threatening.

- Do not climb into the appliance.
- Do not carry out cleaning work in the chamber alone.

⚠ DANGER**Live parts**

When covers are removed, live parts are exposed and contact with these parts may result in electric shock. Electric shock can have serious health consequences including death.

- Only authorised persons may carry out electrical installation work.
- Before starting work, disconnect the unit from the power supply.
- Ensure that the unit is fully de-energised.
- Secure the unit to prevent it from being switched on again.

⚠ WARNING**Poisonous or explosive vapours and gases**

When loading the unit with an unsuitable load, poisonous or explosive vapours or gases may be produced. This could cause the unit to explode, and persons could be severely injured or poisoned.

- The unit may only be loaded with materials and substances which cannot form any toxic or explosive vapours at the set temperature and which cannot explode, burst or ignite.

⚠ WARNING**Overheating of the appliance when door is open**

Leaving the door open during operation can cause the appliance to overheat or pose a fire hazard.

- Do not leave the door open during operation.

2.3 Requirements to be met by Operating Personnel

The appliance may only be operated and maintained by persons who are of legal age and have been instructed accordingly. It is intended to be operated and maintained by trained personnel employed by the owner.

Repairs may only be performed by qualified electricians. The guidelines in the separate service manual must be observed.

2.4 Responsibility of the Owner

The owner of the unit

- is responsible for the flawless condition of the unit and for operating it in accordance with its intended use;
- is responsible for ensuring that persons who operate or service the unit are qualified to do this, have been instructed accordingly and are familiar with these operating instructions;
- must know the applicable guidelines, requirements and operational safety regulations, and train staff accordingly;
- is responsible for ensuring that unauthorised persons cannot access the unit;
- is responsible for ensuring that the maintenance plan is adhered to and that maintenance work is properly carried out;
- has to ensure that the unit and its surroundings are kept clean and tidy, for example through corresponding instructions and inspections;

- is responsible for ensuring that personal protective clothing is worn by operating personnel, e.g. work clothes, safety shoes and protective gloves.

2.5 Product Use

2.5.1 Intended Use

The ULT series of ULT freezers are designed for professional use for the safe storage of temperature-sensitive samples, materials and other items suitable for ultra-low temperature storage within a temperature range of -50 °C to -86 °C. They are used for storage at a defined and constant ultra-low temperature, particularly over extended periods.

The device is particularly suitable for use in laboratories and in the life sciences, research and pharmaceutical sectors. Furthermore, it can be used for similar technical applications where a stable ultra-low temperature is required for samples or materials.

Intended use includes operating the equipment exclusively within the limits specified in the technical data and in these operating instructions. Only suitable, non-hazardous goods that do not damage the equipment and do not pose a risk of fire, explosion, corrosion or any other hazard may be stored.

This device is not a medical device.

2.5.2 Improper Use

Any use of the appliance that deviates from the intended use for ultra-low temperature storage as described in this user manual is considered improper. This applies in particular to operation outside the specified technical data, environmental conditions and safety requirements set out in this operating manual.

In particular, the following constitutes improper use:

- the storage of substances, preparations or materials that are flammable, explosive, corrosive or capable of producing hazardous gases, or that may damage parts of the equipment;
- the storage of goods without appropriate protective measures, if this could endanger people or contaminate the interior of the unit;
- Installation, commissioning, operation, cleaning, maintenance or repair carried out by personnel who have not received training, are not sufficiently qualified or are not authorised;
- operation with a damaged device, a damaged power cable, an obvious malfunction, or with modified, bypassed or missing safety devices;
- the use of spare parts, accessories or consumables not approved by the manufacturer, as well as unauthorised modifications or alterations to the device;
- operation outside the specified installation and ambient conditions, as well as the ingress of water or cleaning agents into the unit.

Any use other than that for which the equipment is intended may result in personal injury or damage to the equipment or stored goods.

2.6 Changes and Alterations

The device must not be modified or altered without authorization. No parts may be added or installed that are not approved by the manufacturer.

Unauthorized modifications or alterations will invalidate the CE conformity of the device and the device must no longer be operated.

The manufacturer shall not be liable for damage, hazards or injuries resulting from unauthorized modifications or alterations or from failure to comply with the regulations in these instructions.

2.7 Behaviour in case of Malfunctions and Irregularities

i

The unit must only be used in a flawless condition. If you, as the operator, notice irregularities, malfunctions or damage, immediately turn off the unit and inform your line manager.

i

You can find information on troubleshooting in the chapter ▶7 Malfunctions, Warning and Error Messages.


See also

📄 Malfunctions, Warning and Error Messages [▶ 32]

2.8 Switching off the Unit in an Emergency

i

If possible, turn off the appliance at the control panel (see ▶6.7.2 Switch off appliance)

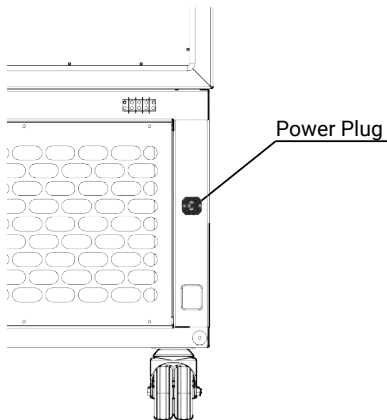- ▶ Unplug the power cord from the power source.
 - ⇒ This disconnects the device from the power supply on all poles.
 - ⇒ The device behaves as it would during a power outage (see ▶7.1.6 Power Failure)

3. Construction and Description

3.1 Design

1	Pressure equalisation	2	Nameplate
3	USB port	4	Castors
5	External door with control panel	6	Interior door
7	Door handle with door lock	8	Refrigeration unit

3.2 Description of Function

The unit can cool the interior to a temperature range between -50 °C and -86 °C. Cooling is provided by a two-stage cascade system. Both stages are fitted with a capillary injection system.

The ULTeco’s compressors are also fitted with an inverter, which regulates the speed – and thus the temperature – in an energy-efficient manner tailored to demand.

3.3 Materials

Components	Material
Housing (door)	Stainless steel 1.4016 – ASTM 430

Components	Material
Housing (top panel + side panels + rear panel)	Stainless steel 1.4301 – ASTM 304
Chamber (incl. cover)	Stainless steel 1.4301 – ASTM 304
Sliding compartments	Aluminium
Door seal	Silicone
Insulation	Plastic insulation boards
ULT battery	Type: ES 12-12 Chemical system: Lead-acid battery Weight: 3.4 kg Capacity: 12 V Labelling: "maintenance-free"
ULTeco battery	Type: LFP 10-12 Chemical system: lithium battery Weight: 1.06 kg Capacity: 12.8 V Labelling: "maintenance-free"

The chamber load of the appliance must be carefully checked for chemical compatibility with the indicated materials.

3.4 Connections and Interfaces

3.4.1 Electrical Connection


This unit is designed for operation on an electrical power system with a maximum system impedance $Z_{\max}$ at the point of transfer (service line) of 0.292 Ohm. The operator must ensure that the unit is only operated on an electrical power system that meets these requirements.

If necessary, ask your local utility company what the system impedance is. Observe the country-specific regulations when making connections (e.g. in Germany DIN VDE 0100 with earth leakage circuit breaker).

3.4.2 Communication Interfaces

The communication interfaces are intended for devices that comply with the requirements of IEC 60950-1.

Ethernet interface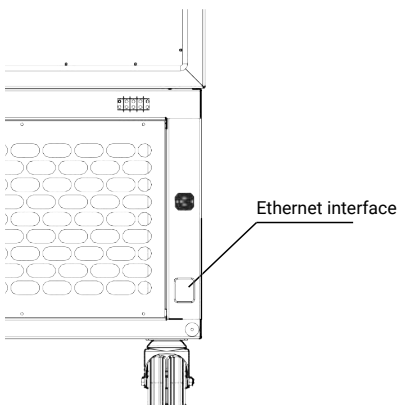

The device can be connected to a network via an Ethernet interface, and logs can be retrieved.

For identification purposes, each unit connected must have its own unique IP address. A description of how to set the IP address is provided in chapter IP Address and Subnet Mask.

The unit can be directly connected to a computer / laptop using an optional USB to Ethernet converter (see ▶3.9 Scope of Delivery).

USB interface

The unit comes with a USB port as standard in accordance with the USB specification. With this you can:

- Logs can be exported from the device to a USB storage device (see Log)
- The USB port is located on the front left-hand side of the device.


See also

 Scope of Delivery [► 17]


3.4.3 Potential-free contact

The device has a potential-free contact to which an external warning system can be connected.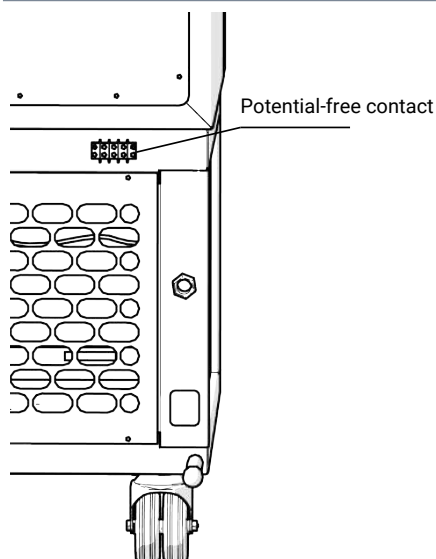

The potential-free contact is located on the right side of the back of the device.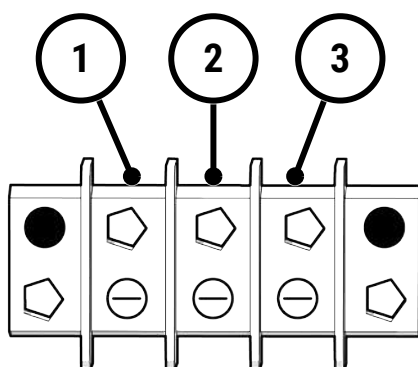

1 C = Common

2 NC = Normally closed

3 NO = Normally open

- Contact C must be connected to the power supply.
- By default, current flows between contact C and contact NC when contact NC is closed.
- With contact NO, no current flows by default.

The following fault messages are displayed via the dry contact (see ►7.1.2 Fault Messages):

- E1
- E2
- E4
- LO1
- HI1
- HI2

- Simulate the alarm on the cabinet to identify the correct contact (NO or NC).

3.5 Nameplate

memmert GmbH+Co.KG

Äußere Rittersbacher Str. 38
 D91126 Schwabach, Germany
 Made in the Netherlands

3 SYSTEM

4 MODEL

5 SERIAL NO.

6 TYPE REFRIGERANT:

7 CAPACITY

8 VOLTAGE

Imax

: cascade cooling

: ULT360eco

: U336.0001

Stage 1 : R290 = 310 gram

Stage 2 : R170 = 136 gram

: 360 liter

: 230 V / 50-60 Hz. - 1 ph



: 3,92 A

1	Contact details	2	Test marks and warning symbols
3	Refrigeration system	4	Model name
5	Serial Number	6	Refrigerant
7	Interior volume	8	Electrical data

3.6 Technical Data

Unit size				240	360	500	750	1000
Stainless steel interior	Volume		l	244	360	506	750	966
	Width	A	mm	450	450	600	890	1035
	Height	B	mm	1111	1276	1345	1345	1276
	Depth	C	mm	488.5	626.5	626.5	626.5	731.5
	Max. loading per shelf		kg	140	160	180	200	220
	Max. loading per unit		kg	280	320	360	400	440
Patterned stainless steel housing	Width	D	mm	700	700	850	1140	1285
	Height	E	mm	1832	1992	1999.5	1999.5	1992
	Depth	F	mm	738	888	888	888	973
Temperature	Setting temperature range		°C	-50 up to -86				
	Adjustment precision		°C	0.1				
Further data	Power consumption (50/60 Hz) ULT	230 V	W	589	684	960	1130	1375
	Max. current consumption (50/60 Hz) ULT	230 V	A	3.5	3.6	4.8	5.4	7.0
	Power consumption (50/60 Hz) ULTeco	230 V	W	470	500	540	560	600

Unit size				240	360	500	750	1000
	Max. current consumption (50/60 Hz) ULTeco	230 V	A	3.35	3.92	4	4.21	4.7
	Noise level ULT		dBA	51	52	53	54	55
	Noise level ULTeco		dBA	49	49	49	49	49
Packaging data	Net weight		kg	245	285	343	380	425
	Gross weight		kg	310	370	430	550	590
	Width		mm	900	1080	1080	1410	1440
	Height		mm	2050	2220	2220	2220	2250
	Depth		mm	900	1080	1080	1080	1080

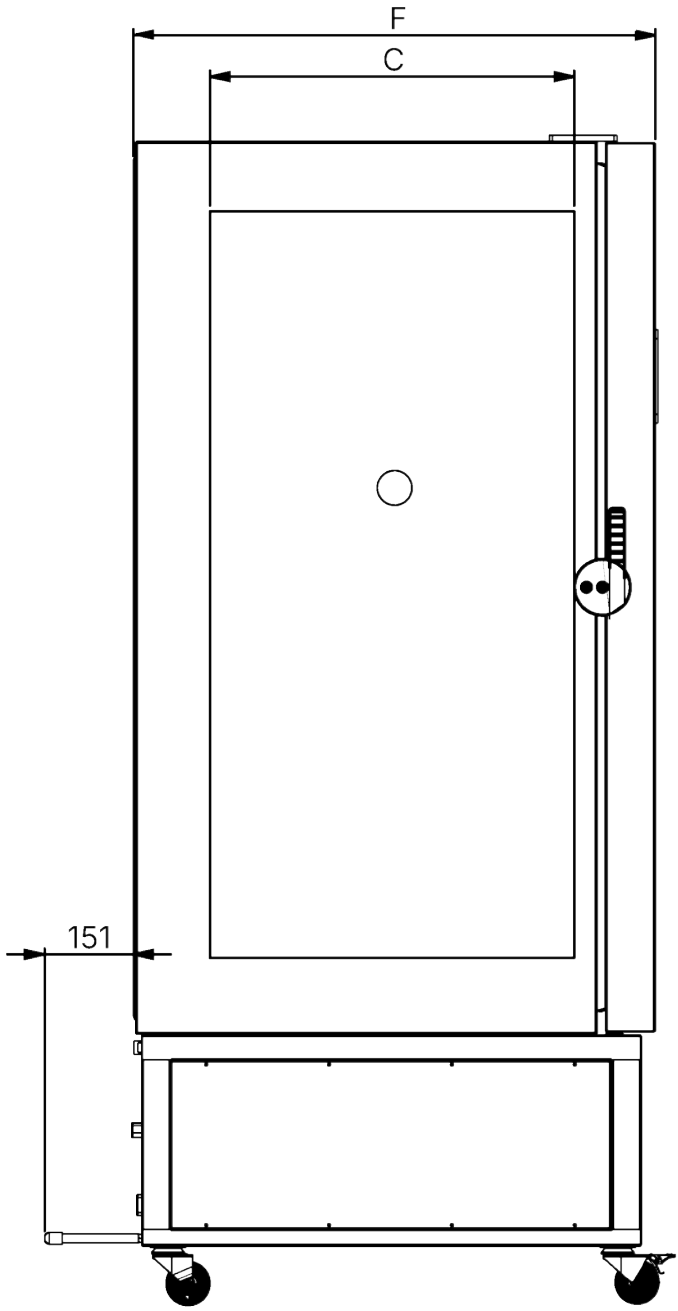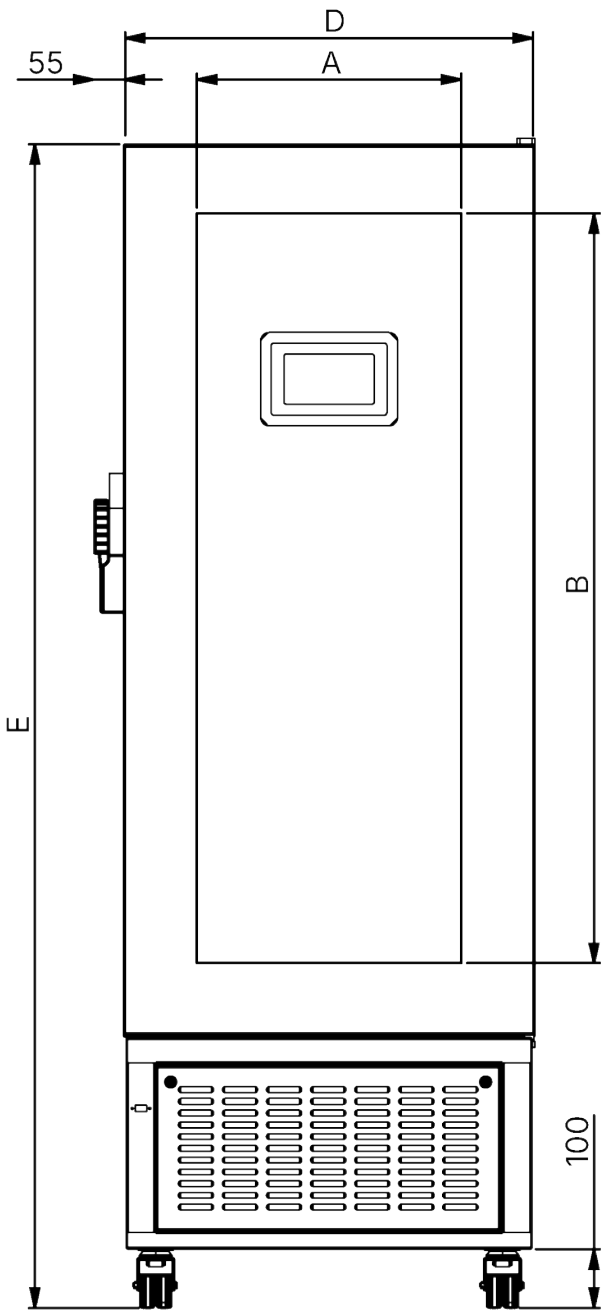Information on the refrigerant ULT

Information on the refrigerant ULT		240	360	500	750	1000
First-stage refrigerant		R455A	R455A	R455A	R455A	R455A
Second-stage refrigerant		R170	R170	R170	R170	R170

Information on the refrigerant ULT		240	360	500	750	1000
Fluid group		1	1	1	1	1
Filling capacity, first stage	g	450	450	650	700	750
Filling capacity of the second stage	g	150	160	180	190	230
Cooling capacity	W	589	684	960	1130	1375
Max. ambient temperature	°C	28	28	28	28	28
Max. operating pressure	bar	22.5	22.5	22.5	22.5	22.5
Safety class		A3	A3	A3	A3	A3

Information on the ULTeco refrigerant

Information on the ULTeco refrigerant		240	360	500	750	1000
First-stage refrigerant		R290	R290	R290	R290	R290
Second-stage refrigerant		R170	R170	R170	R170	R170
Fluid group		1	1	1	1	1
Filling capacity, first stage	g	310	310	330	360	360
Filling capacity of the second stage	g	130	136	146	160	170
Cooling capacity	W	470	500	540	560	600
Max. ambient temperature	°C	28	28	28	28	28
Max. operating pressure	bar	22.5	22.5	22.5	22.5	22.5
Safety class		A3	A3	A3	A3	A3

3.7 Applied Directives and Standards

Based on the standards and guidelines listed below, the products described in this manual carry a CE mark from Memmert:

- Directive 2014/30/EU with amendments (Council Directive on the approximation of the laws of the Member States relating to electromagnetic compatibility). Standard complied with: EN 60601-1-2, as published in the Official Journal of the European Union.
- Directive 2014/35/EU (Directive of the European Parliament and of the Council of 26 February 2014 on the harmonisation of the laws of the Member States relating to the making available on the market of electrical equipment designed for use within certain voltage limits). Standard complied with: EN 61010-1, as published in the Official Journal of the European Union.

3.7.1 Material Compliance

We confirm that we always draw the attention of our suppliers to the legal restrictions on materials in accordance with our **Company Standard for Material Compliance of Memmert GmbH + Co KG** to ensure they take the original publications by the legislative authority into consideration at all times. The suppliers and deliveries must comply with all material compliance requirements which are relevant or specified in the company standard. By taking this approach, and by making our own observations, we are always able to stay abreast of developments to the best of our knowledge and ability.

The latest material compliance statement can be accessed via the link <https://www.memmert.com/de/downloads>.

3.8 Ambient Conditions

- The unit must only be used in closed rooms and in the ambient conditions listed below:

Ambient temperature	+15 °C to +28 °C
Air humidity	max. 70% non-condensing
Overvoltage category	II
Contamination level	2
Installation altitude a.s.l.	max. 2000 m a.s.l.

- The unit may not be used in Ex zones. The ambient air must not contain explosive dusts, gases, vapours or gas-air mixtures. The unit is not explosion-proof.
- Heavy dust production or aggressive vapours in the vicinity of the unit could lead to sedimentation in the interior and, as a consequence, could result in short circuits or damage to electrical parts. For this reason, sufficient measures to prevent large accumulations of dust or aggressive vapours should be taken.

3.9 Scope of Delivery

- Mains connection cable
- 2 spacers
- Operating manual
- 3 drawers
- Key for door lock

4. Delivery, Transport and Setting Up

4.1 Safety

⚠ CAUTION

Lifting the appliance incorrectly

The appliance is heavy. The appliance is heavy, so you could injure yourself if you try to lift it on your own.

- Only move the unit using a pallet truck or forklift.

240	360	500	750	1000
				

⚠ CAUTION

Crushing hazard due to heavy equipment

The appliance is heavy. Crushing injuries to hands or feet can occur when transporting and installing the unit.

- Wear protective gloves and safety boots.

⚠ CAUTION

Risk of injury due to the appliance falling over during transport

The appliance is heavy. The appliance could fall over and seriously injure you.

- Never tilt the appliance and only transport it in the upright position without load (except for standard accessories such as grids or shelves).
- Appliances with castors always have to be moved by at least two people.

4.2 Delivery

The appliance is supplied packed in cardboard on a wooden pallet.

4.3 Transport

The unit can be transported in different ways depending on its size:

- With a forklift truck or pallet truck; move the forks of the truck entirely under the pallet
- If the appliance has its own castors, also release the locking device on the (front) castors

4.4 Unpacking

- Do not unpack the appliance until you reach the installation site.
- Remove the cardboard packaging by pulling it upwards or carefully cutting along an edge or unscrew and remove wooden crate.

Checking for completeness and transport damage

- Check the delivery note to ensure the delivery is complete.
- Check the unit for damage.

If you notice deviations from the scope of delivery, damage or anything unusual, do not put the unit into operation and inform the haulage company and the manufacturer.

Remove the protective film from the housing

- Remove the protective film from all components before commissioning.

Remove transport securing devices

- Remove the transportation lock. It is located between the door hinge, door and frame and has to be removed after opening the door.

Recycling of packaging material

- Dispose of the packaging material (cardboard, wood, foil) in accordance with the applicable disposal regulations for the respective material in your country.

4.5 Storage after Delivery


If the unit is to be stored initially after delivery:

- Observe the storage conditions (see ►9 Storage and transport).


4.6 Setting Up

4.6.1 Preconditions

- ✓ The distance between the wall and the rear panel of the appliance must be at least 15.5 cm.
- ✓ The clearance from the ceiling must not be less than 15 cm and the side clearance from walls or nearby appliances must not be less than 15 cm. Sufficient air circulation in the vicinity of the appliance must be guaranteed at all times.
- ✓ For appliances with castors or for appliances standing on sub frames with castors, always position the castors in a forward direction and lock the castor brakes to ensure that the appliances remain securely in place.
- Place the unit in the designated position as shown below.

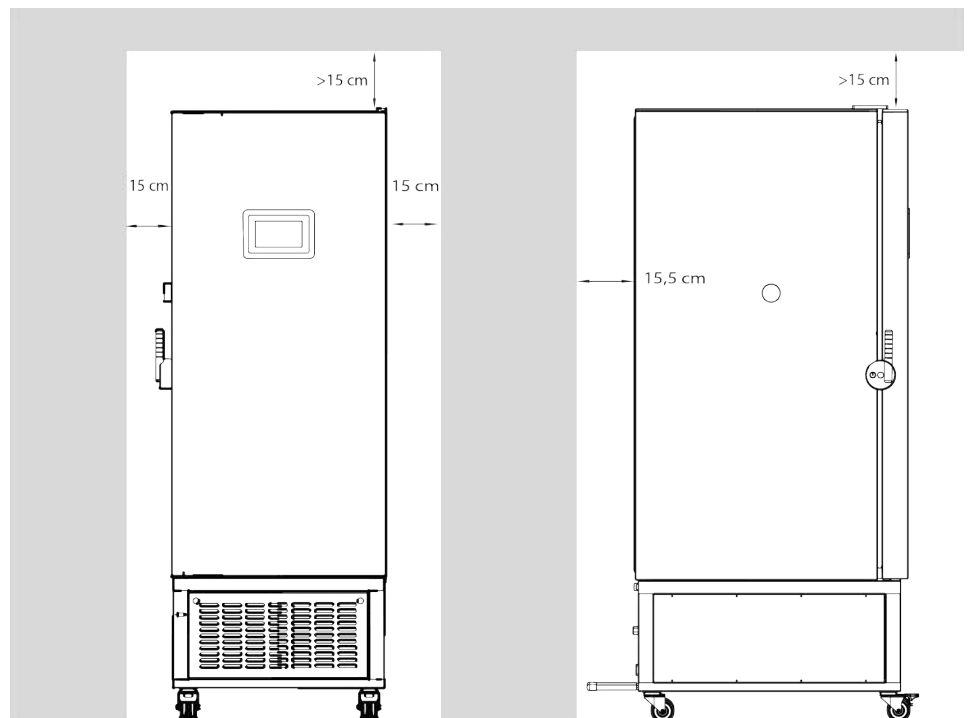
4.6.2 Spacer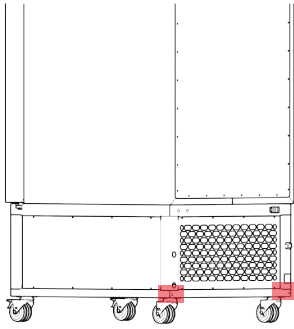

- The spacers are not pre-assembled at the factory.
- The spacers must be fitted to the rear panel before commissioning.

-
1. Screw the spacers into the threads on the back of the unit.

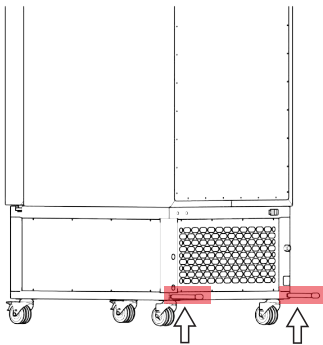5. Putting into Operation

5.1 Putting into Operation for the First Time

⚠ WARNING

Condensation in the electrical components may cause short circuits.

Due to temperature fluctuations during transport, condensation may form inside the unit.

- After transporting or storing the unit in humid conditions, remove it from its packaging and allow it to acclimatise for at least 24 hours in normal ambient conditions.
- Do not connect the unit to the power supply during this time.

NOTICE

When putting the unit into operation for the first time, do not leave it unattended until it has reached a steady state.

- Please observe the national regulations when connecting the unit.
- Observe the connected loads and power ratings (see ▶3.5 Nameplate and ▶3.6 Technical Data).
- Be sure to establish a safe PE conductor connection.

See also

- 📄 Nameplate [▶ 14]
- 📄 Technical Data [▶ 14]


5.2 Connecting the Unit to the Power Supply

Observe the country-specific regulations when making connections (e.g. DIN VDE 0100 with earth leakage circuit breaker, in Germany).

Observe the connected loads and power ratings (see ▶3.5 Nameplate and ▶3.6 Technical Data).

Be sure to establish a safe PE conductor connection.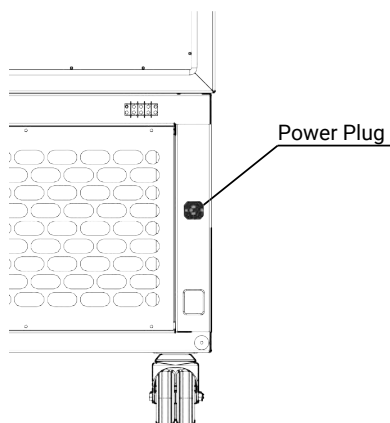


Route the power cable so that

- nobody can trip over it.
- it cannot come into contact with any hot parts.
- it is easily accessible at all times and the plug can be pulled out quickly in the event of a fault or emergency, for example.

See also

- 📄 Nameplate [▶ 14]
- 📄 Technical Data [▶ 14]

5.3 Switching on Unit

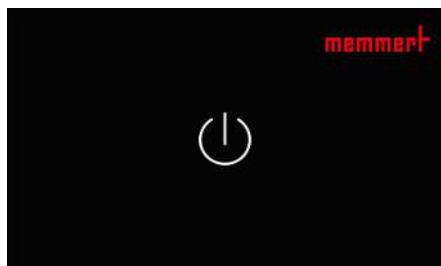- ✓ The device must be connected to the electricity grid.
 - 1. Tap the power icon to switch on the device.
-

6. Operation and Control

⚠ CAUTION

Slip hazard

Small amounts of condensed water may leak from the appliance when it is in operation.

- Always wear shoes with non-slip soles and wipe up the condensed water immediately.

⚠ CAUTION

Injuries from cold surfaces

The surfaces of the chamber and chamber load may be very cold.

- Always wear gloves when you reach into the appliance.

6.1 Operating Personnel


The appliance may only be operated by persons who are of legal age and have been instructed accordingly. Personnel who are to be trained, instructed or who are undergoing general training may only work with the appliance under constant supervision of an experienced person.

6.2 Opening the Door

1. Open the door by pulling the door handle downwards.

6.3 Switching on Unit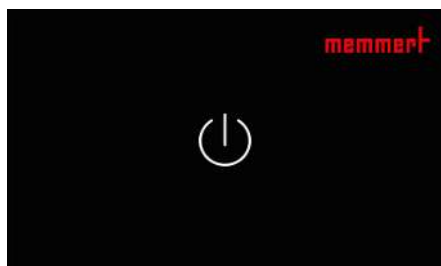


- ✓ The device must be connected to the electricity grid.

1. Tap the power icon to switch on the device.

6.4 Role-based access control system

Role	Adjustable parameters
Administrator	all
User	30, 31, 80, 81, 82, 83, 84, 90, 91, 95, 99

6.5 Log in

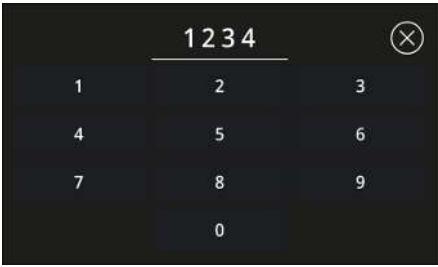
Once the device has been switched on, the login window will appear.

- 1. Enter your four-digit login code.
⇒ You will be logged in automatically.

- The factory default login code for the **User** role is 1234.
- The factory-set login code for the **Administrator** role is 5214.
- The factory reset code is 6525.

6.6 Dashboard

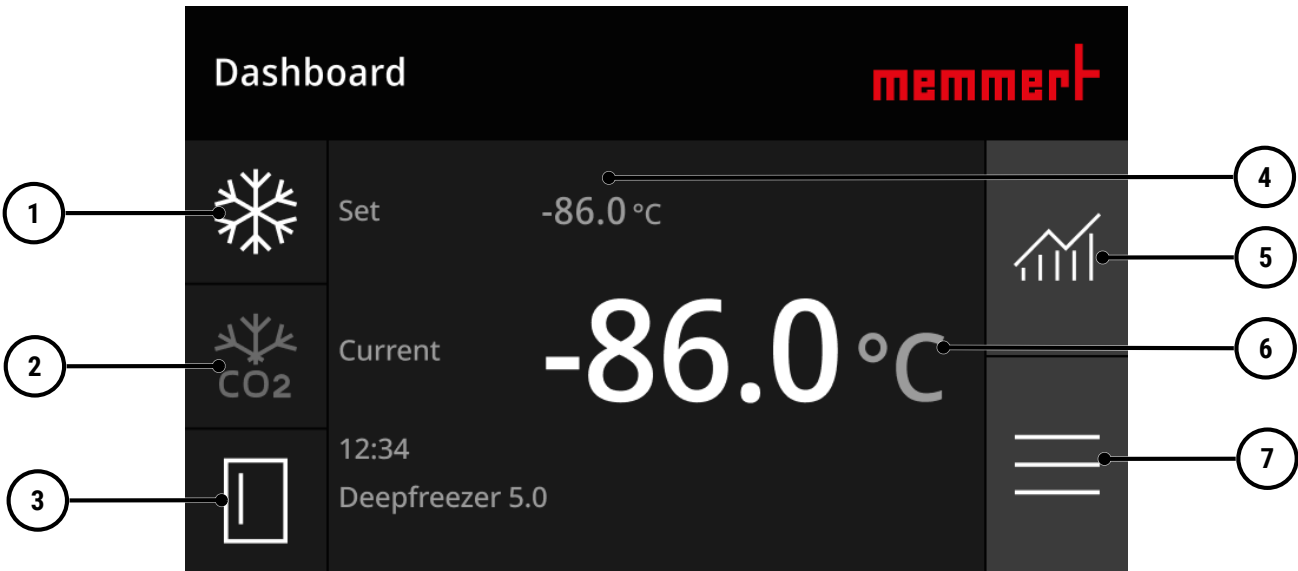1	Adjusting temperature	2	CO ₂ emergency cooling indicator (optional)
3	Door open indicator	4	Setpoint temperature
5	Display temperature graph	6	Actual temperature
7	Menu [► 26]		

6.6.1 Distance View

After 4 minutes of inactivity, the control unit switches to Distance View.


1. Tap the control unit to exit Distance View.

After using Distance View, you will need to log in again.


2. Select the padlock icon to log in.

6.6.2 Adjusting temperature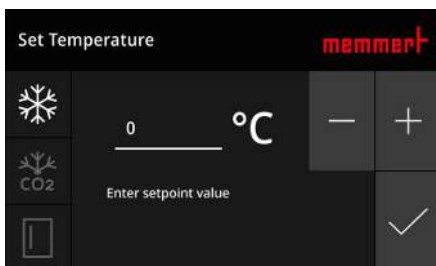


1. On the dashboard, select **Set temperature**.
2. Use the plus and minus symbols to set the desired target temperature.
3. Select the tick icon to confirm
 - ⇒ The appliance begins the cooling process.

6.6.3 Display temperature graph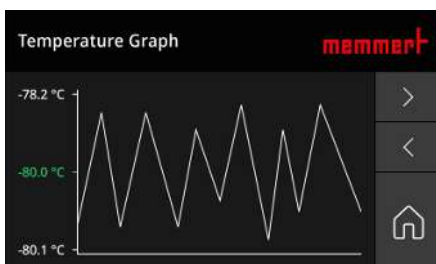

- Use the arrows to view the temperature graph for different time periods.
- The house icon takes you back to the dashboard.

6.7 Menu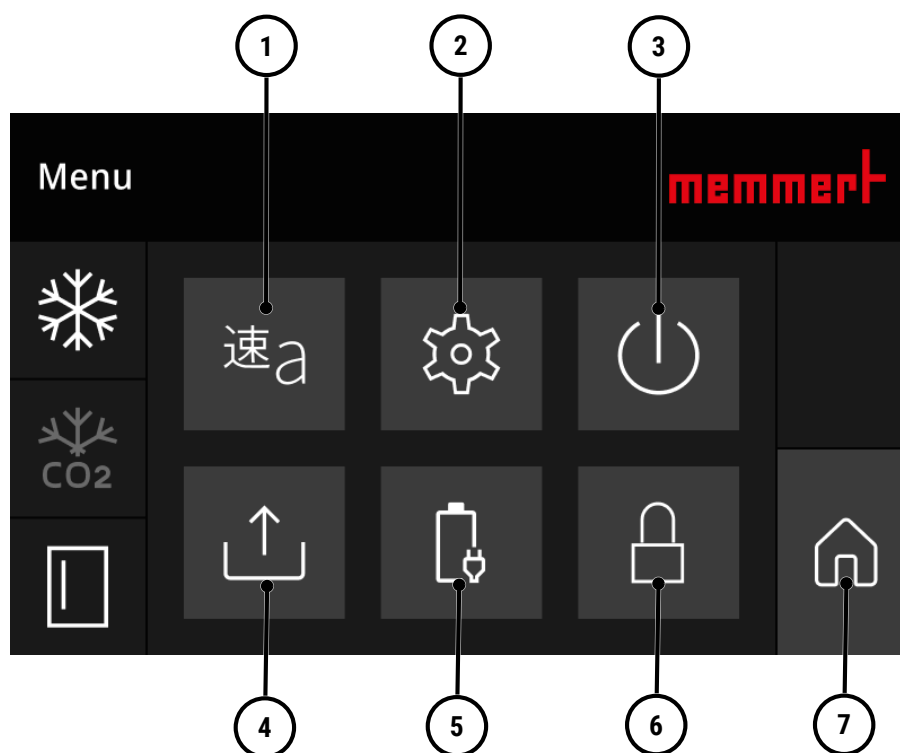

1	Language settings	2	Options [► 28]
3	Switch off appliance	4	Transfer log data
5	View battery information	6	Lock device
7	Back to the dashboard		

6.7.1 Language settings


The languages available are English, German, Italian and Dutch.

1. In the menu, select **Language Settings**.
2. Select the desired interface language.
⇒ The selected language is highlighted in dark.
3. Tap the three-line icon to return to the menu.

6.7.2 Switch off appliance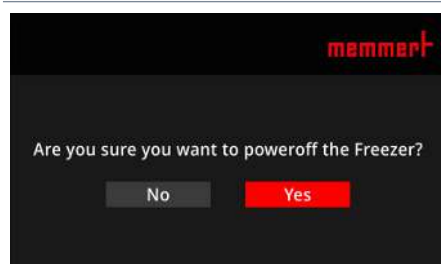

1. From the menu, select **Turn Off Device**.
2. Confirm in the dialog box.
⇒ The device turns off.

6.7.3 Transfer log data


The unit continually logs all relevant measured values, settings and error messages at 1-minute intervals. The internal log memory is a continuous memory type. The log function cannot be switched off and is always active. The measurement data is stored in the unit, safe from manipulation. If the power supply is interrupted, the time of the power failure and when the power was restored are stored in the unit.

Log data from various time periods can be downloaded to a USB storage device via the USB interface or via Ethernet.


The log memory of the appliance is not modified or deleted by reading it out.

- The USB port is located on the front left-hand side of the device.
1. Insert the USB storage device into the USB port.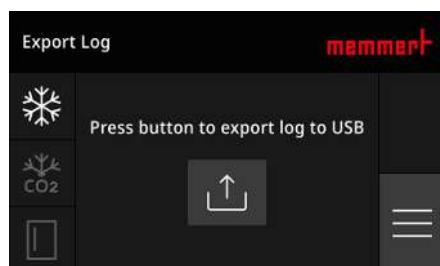


2. Select **Transfer log data** from the menu.
3. Select the export icon to transfer the log data.
 - ⇒ The log data is transferred to the USB storage device.
4. Tap the three-line icon to return to the menu.

6.7.4 View battery information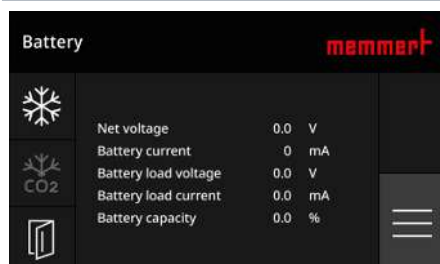


- This screen displays the current battery status information.
- The three-bar icon takes you back to the menu.

6.8 Options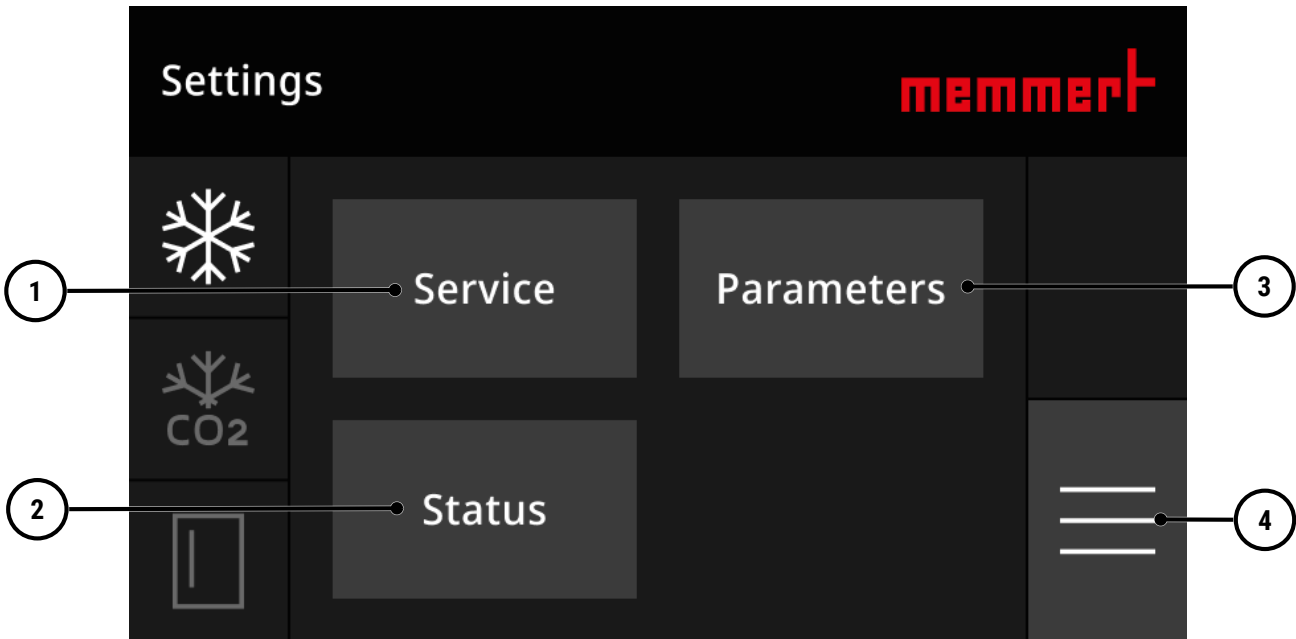


1	Record service activities	2	Display temperature and compressor status
3	Setting Parameters	4	Back to the menu

6.8.1 Record service activities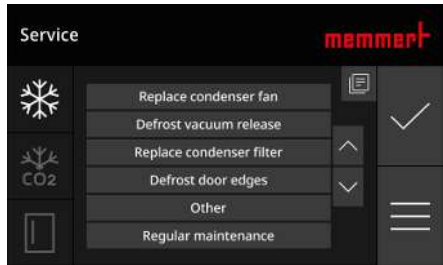


- This is where details of the service work carried out can be recorded.
1. In the settings, select **Log service activities**.
 2. Use the arrow keys to select the service task that has been carried out.
 3. Select the tick icon to confirm.
⇒ The service work carried out is recorded in the log.
 4. Tap the three-line icon to return to Settings.

6.8.2 Display temperature and compressor status

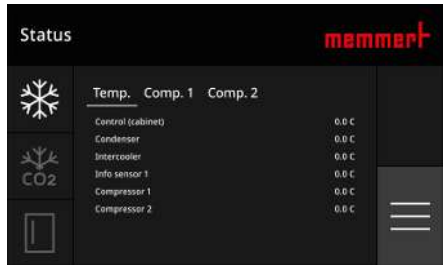
- Temperature status information is displayed here.
- The three-bar icon takes you back to the settings.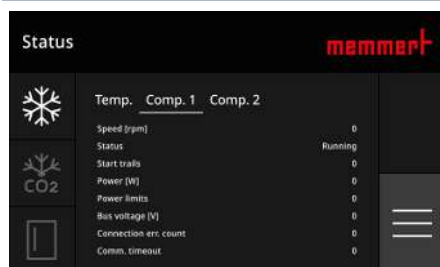

- This is where the compressor status information is displayed.
- The three-bar icon takes you back to the settings.

6.8.3 Setting Parameters

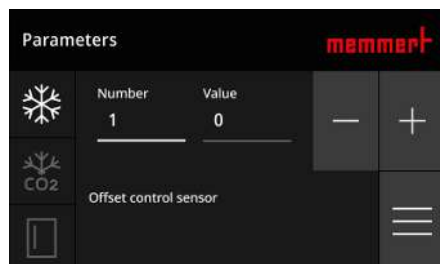1. Select the parameter number using the plus and minus symbols.
2. Set the parameter value using the plus and minus symbols.
3. Tap the three-line icon to return to Settings.

Parameter

No.	Definition	Role	Setting range	Unit	Ex works
Group 0					
1	Offset sensor	Administrator	-150.0 .. 150.0	K	2.0
2	Offset sensor for the condenser	Administrator	-150.0 .. 150.0	K	0.0
3*	Capacitor probe display	Administrator	0.0 .. 0.0	°C	0.0
4	Freezer type (0=upright freezer, 1=chest freezer)	Administrator	0 .. 1	-	0
Group 1					
10	Minimum setpoint	Administrator	-100.0 .. 25.0	°C	-86.0
11	Maximum setpoint	Administrator	-100.0 .. 25.0	°C	-50.0
12	Display (0 = for 0.5 °C, 1 = for 1.0 °C)	Administrator	0 .. 1	-	0
Group 2					
20	Temperature difference	Administrator	-150.0 .. -0.1	°C	-0.6
21	Start-up delay following a power cut	Administrator	0 .. 99	Min	0
22	Delay in the fan switching off	Administrator	0 .. 9999	Sec	300
Group 3					
30	Minimum alarm temperature (control sensor)	User	-100.0 .. 25.0	°C	-85.0
31	Maximum alarm temperature (sensor)	User	-100.0 .. 25.0	°C	-70.0
32	Maximum alarm temperature (condensation sensor)	Administrator	0.0 .. 100.0	°C	38.0
33	Activate the alarm relay when the maximum alarm is triggered	Administrator	0 .. 1	-	1
34	Delay of the door-opening alarm	Administrator	0 .. 999	sec.	30
37	Alarm temperature relay	Administrator	-100.0 .. 25.0	°C	-45.0
38	Minimum alarm delay (Lo)	Administrator	0 .. 999	Sec	0
39	Maximum alarm threshold (Hi)	Administrator	0 .. 999	Sec	0
Group 4					
40	Battery capacity	Administrator	1 .. 150	Ah	10
46	Start delay for compressor 2	Administrator	0 .. 1800	Sec	60
Group 5					
50	Start-up temperature for CO ₂ cooling	User	-100.0 .. 25.0	°C	-60.0

51	Pulse CO ₂ emergency cooling	Administrator	1 .. 999	Sec	5
52	Pause in emergency CO ₂ cooling	Administrator	1 .. 999	Sec	20
53	CO ₂ emergency cooling enabled when mains power is available (0=Disabled, 1=Enabled)	Administrator	0 .. 1	-	1
Group 6					
60	Slow temperature difference	Administrator	0.0 .. 25.0	°C	6.0
61	Speed Temperature Slow	Administrator	1 .. 100	-	20
62	Time until the next service	Administrator	-9999 .. 9999	-	365
63	Logo	Administrator	0 .. 6	-	0
64	Setpoint for service delay	Administrator	0 .. 9999	-	365
Group 7					
70	Alarm relay control	User	0 .. 999	Sec	60
78	Battery voltage monitoring off	Administrator	0 .. 1	-	0
79	Battery voltage too low	Administrator	0.0 .. 999.9	volt	11.5
Group 8					
80	Current date: Year	User	0 .. 9999	-	0
81	Current date: Month	User	1 .. 12	-	0
82	Current date: Day	User	1 .. 31	-	0
83	Current time: hour	User	0 .. 23	-	0
84	Current time: minutes	User	0 .. 59	-	0
Group 9					
90	Access code	User	1111 .. 5555	-	1234
91*	Network address (IP)	User	0 .. 0	-	1
95*	SD card software version	User	0.00 .. 0.00	-	0.00
96*	Serial Number	Administrator	0 .. 0	-	0
97*	Date created (YY-DD) (YY is the year, DD is the day of the year, with 1 representing 1 January)	Administrator	0 .. 0	-	0
99	Restore factory settings (2119) (Settings & IP) (automatic restart)	User	1111 .. 9999	-	0
Group 20					
203	Offset of the charge air sensor	Administrator	-50.0 .. 10.0	K	0.0
204	Disable UPS main voltage fault (0=no, 1=yes, 2=no automatic reset)	Administrator	0 .. 2	-	0
205	Disable compressor error (0=no, 1=yes)	Administrator	0 .. 1	-	0
206	Disable Summer (0=no, 1=yes, 2=no, including E25)	Administrator	0 .. 2	-	0
207	Reset all alarms automatically (0=no, 1=yes)	Administrator	0 .. 1	-	0
Group 21					
210	Max. compression speed 1	Administrator	0 .. 5000	-	5000
211	Max. speed of compressor 2	Administrator	0 .. 5000	-	5000
213	Activate Stage 2 intercooler temperature	Administrator	-999.9 .. 999.9	°C	-28.0
214	Deactivation of intercooler temperature stage 2	Administrator	-999.9 .. 999.9	°C	-15.0
215	E1 Compressor 1 speed (1800 rpm .. 210 rpm)	Administrator	0 .. 100	%	0
216	E1 Variable-speed compressor 2 (1800 rpm .. P211 rpm)	Administrator	0 .. 100	%	0
217	Minimum compressor speed	Administrator	0 .. 4500		1800
Group 22					
220	Trap 1 P-band	Administrator	0.0 .. 25.0	°C	5.0
221	Scenario 1: I-time	Administrator	0 .. 150	Min	5

222	Trap 2 P-Band	Administrator	0.0 .. 25.0	°C	5.0
223	Case 2: I-time	Administrator	0 .. 150	Min	15
224	Case 2 D-factor / 100	Administrator	0 .. 100	-	0
Group 23					
231	Cool-Step output difference	Administrator	0 .. 100	%	3
232	Compr. 2 Allow deactivation if the SP is above this value	Administrator	-999.9 .. 999.9	°C	-75.0
233	Intercooler setpoint	Administrator	-999.9 .. 999.9	°C	-32.0
234	Restart Initial phase Setpoint	Administrator	-999.9 .. 999.9	°C	-50.0
Group 24					
240	Setpoint temperature	Administrator	-100.0 .. 25.0	°C	0.0
242	Start status	Administrator	0 .. 9999	-	0
Group 25					
250	Analogue output for cabinet temperature at 0 V or 0 mA	Administrator	-999.9 .. 999.9	°C	0.0
251	Analogue output for cabinet temperature at 10 V or 20 mA	Administrator	-999.9 .. 999.9	°C	-100.0
Group 26					
260	Apply the cooling difference (P20) to stop the inverter compressor	Administrator	0 .. 1	-	1
261	Delay in restarting the compressor after it has been shut down by P260	Administrator	0 .. 9999	sec.	10
Group 99					
992	Restart the controller (1111)	Administrator	1111 .. 9999	-	0
993	Freezer on/off (0=off, 1=on)	Administrator	0 .. 1	-	0
995	Language (0=EN, 1=NL, 2=DE, 3=IT)	Administrator	0 .. 99	-	0
996*	Software version of the UPS board	User	0.00 .. 0.00	-	0.00
997	Software version of the UPS board	User	0.00 .. 0.00	-	0.00

Parameters marked with an * are read-only.

Note: Some parameters are only available for ULT eco.

6.9 Defrost the unit

1. Turn off the unit (see ►6.7.2 Switch off appliance).
2. Open the door (see ►6.2 Opening the Door).
3. Remove all shelves.
4. Leave the door open.
5. If necessary, place towels or similar items underneath to catch any dripping liquid.
6. Wait until the appliance is completely free of ice.

7. Malfunctions, Warning and Error Messages

⚠ DANGER

Risk of electric shock from unauthorised troubleshooting

Errors requiring intervention inside the unit may only be rectified by qualified electricians.

- Follow the measures listed in the event of a malfunction.
- Contact Memmert International After Sales.

Do not try to rectify appliance errors yourself; instead you should contact Memmert International After Sales or an authorised customer service point.

In case of enquiries, please always state the model and appliance number on the nameplate (see ▶3.5 Nameplate).

See also

- ▢ Nameplate [▶ 14]

7.1 Fault Messages

7.1.1 Dealing with fault reports

- The fault message is also indicated by an audible alarm.

1. Tap the bar displaying the fault message to temporarily disable the beep until the next fault occurs

On the control panel, the fault message is displayed as a red bar at the top.


In Distance View, the fault message is displayed as a red indicator.

Acknowledge fault reports


- ✓ You are logged in.

1. Tap the error message, which is displayed as a red banner at the top of the screen.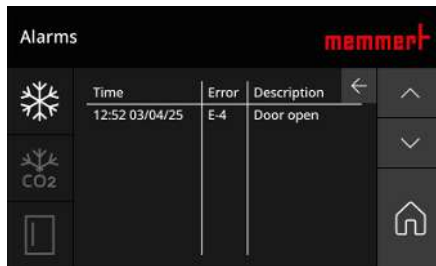

2. Use the arrows in the fault menu to select the fault message you wish to acknowledge.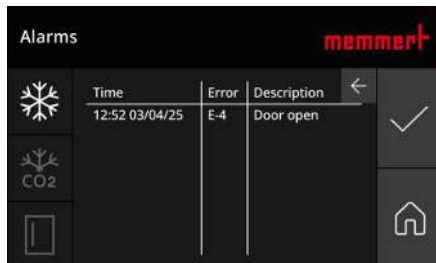


3. Tap the tick to confirm.

⇒ The fault report has been acknowledged.

7.1.2 Fault Messages

Fault report	Cause of the fault	Troubleshooting
E	Faulty condenser sensor	Check the connections
E1	Faulty interior temperature sensor	Check the connections
E2	Mains power is missing	- Check whether there is a power cut - Check the internal fuse
E3	Battery not connected	Check connections, fuses and the battery
E4	No power supply	Check connections and fuses
E5	Battery voltage too low	Check the battery
E6	Door open notification	Close door
E10	Fluctuations in the mains voltage	Notify customer service
E20	Emergency mode active	Notify customer service
Lo1	The temperature inside the vehicle is too low	Notify customer service
Hi1	The temperature inside is too high	Notify customer service
Hi1	Temperature too high after the condenser	Notify customer service

7.1.3 Door open notification

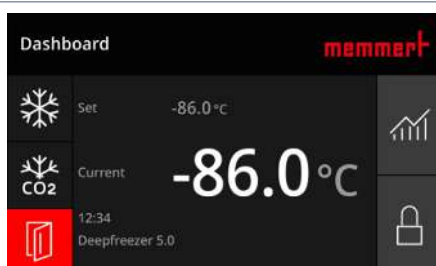The appliance is fitted with a door-open sensor. The 'Door Open' message is displayed in red in the menu.

7.1.4 Emergency mode

If there are any problems or errors in the software, the device switches to emergency mode and displays error message E20. In emergency mode, the unit continues to cool.

1. Please contact customer services.

7.1.5 Connection error

Could not establish connection to the mainboard.
Please check connection.

If the control unit cannot establish a connection to the power unit, an error message will be displayed in full screen.

1. Please contact customer services.

7.1.6 Power Failure

In the event of a power cut, the device will behave as follows:

- The device displays error code E2.
The device runs on battery power for up to 48 hours
- to operate the control unit,
- to continue to log all relevant measurement values, settings and error messages,
- if available, to activate the CO₂ emergency cooling system.

8. Maintenance and Servicing

DANGER

Danger of suffocation inside the appliance

If the appliances is a certain size, you can get accidentally locked in, which is potentially life-threatening.

- Do not climb into the appliance.
- Do not carry out cleaning work in the chamber alone.

DANGER

Live parts

When covers are removed, live parts are exposed and contact with these parts may result in electric shock. Electric shock can have serious health consequences including death.

- Only authorised persons may carry out electrical installation work.
- Before starting work, disconnect the unit from the power supply.
- Ensure that the unit is fully de-energised.
- Secure the unit to prevent it from being switched on again.

CAUTION

Danger of cuts due to sharp edges

Touching sharp edges on the unit may result in cuts.

- Wear protective gloves during all work.
- Be careful when handling sheet metal parts.

8.1 Cleaning

Interior and metal surfaces

Regular cleaning of the easy-to-clean interior prevents build up of material residues that could impair the appearance and functionality of the stainless steel chamber over time.

The metal surfaces of the appliance can be cleaned with normal stainless steel cleaning agents. Make sure that no rusty objects come into contact with the interior or with the stainless steel housing. Rust deposits can lead to an infection of the stainless steel. If rust spots appear on the surface of the interior due to impurities, the affected area must be immediately cleaned and polished.

Plastic parts

Do not clean plastic parts of the appliance with caustic or solvent-based cleaning agents.


Cooling unit

In order to guarantee proper functioning and a long lifetime of the refrigeration unit, it is absolutely essential to remove dust deposits from the cooling unit and the base at least twice a year – the fins and fans must be cleaned with a paintbrush, bottle brush or vacuum cleaner

Observe the specified ambient conditions (see ►3.8 Ambient Conditions) for the installation site of your appliance. If the ambient conditions are unfavourable and there is a lot of dust and dirt, the cleaning intervals should be adjusted as required.

Pull the front panel forwards using both hands.

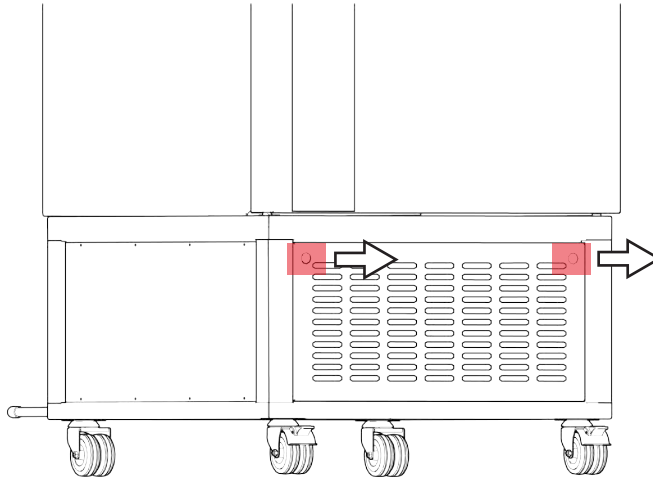**See also**

■ Ambient Conditions [► 16]

8.2 Regular Maintenance

- The appliance must be defrosted at least once a year. However, in conditions of high humidity and frequent opening of the door, defrosting may be required much more often.
 - The filter on the air intake should be checked and cleaned regularly. Cleaning must be carried out every 3 months, or more frequently in dusty environments.
- i**

 - Regular, documented inspections, maintenance and cleaning are essential for the ULT series of freezers to ensure they operate as intended. Failure to carry out and document inspections, maintenance or cleaning may, in certain circumstances, result in the loss of the warranty.

8.3 Repairs and Service

Maintenance and service work may only be carried out by the manufacturer and qualified service providers.

NOTICE

Repairs and service work are described in a separate service manual.

9. Storage, Transport and Disposal

9.1 Storage and Transport

The appliance may only be stored and transported under the following conditions:

- in a dry enclosed, dust-free room
- disconnected from the power supply

9.2 Disposal

This product is subject to Directive 2012/19/EC on Waste Electrical and Electronic Equipment (WEEE) of the European Parliament and EU Council of Ministers. This unit was placed on the market after 13 August 2005 in countries which have already integrated this Directive into their national laws. It must not be disposed of as normal household waste. For disposal, please contact your dealer or the manufacturer. Any units that are infected, infectious or contaminated with materials hazardous to health are excluded from return. Please also observe all other regulations applicable in this context.

Before disposing of the appliance, please render the door locking mechanism unusable, for example to prevent playing children playing with the appliance and being locked inside.

The ULT freezer is fitted with a battery for emergency mode. Remove the battery and dispose of it in accordance with the relevant national regulations.

1. Switch off the appliance.
2. Unplug the mains plug.
3. Pull the front panel forwards using both hands.
4. Disconnect the battery.
5. Lift the battery out upwards.

- Note for Germany:
Do not dispose of the device at public or local collection points

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